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Description:

Test strips for determining the silver concentration and the pH value in fixing baths.

Pack content:

1 aluminum container with 100 test strips

Measuring range:

pH 4–8

Silver: 0–10 g/L Ag⁺

Color gradation:

pH 4 · 5 · 6 · 7 · 8

0 · 0.5 · 1 · 2 · 3 · 5 · 7 · 10 g/L Ag⁺

Additional indications:

1. Ag-Fix cannot be used for fixing baths with pH values lower than 4.
2. To achieve a more exact determination of the silver content in concentrated solutions (more than 5 g silver per liter), dilute the solution with water in a ratio of 1:1 or 1:2. This will move the silver concentration value to the lower range of the test strip, where the reaction is the most sensitive.

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test fields.

Hazard warnings:

This test does not contain hazardous substances that must be labelled.

Instructions for use:

1. Dip the test strip shortly in the fixing bath.
2. Shake off excess liquid.
3. Wait 30 seconds.
4. Compare with the color scale.

Depending on the silver concentration, the yellow test field furthest away turns brown. The pH value can be determined at the same time.

Interferences:

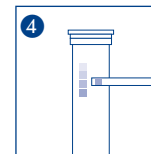
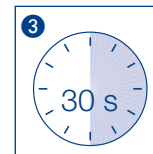
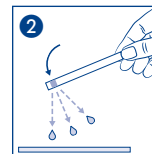
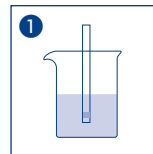
Iron and copper ions cause interference and can cause a false positive reaction. Iron ions lead to these disturbances from a concentration of 1000 mg/L and copper ions from 10 mg/L. These ions can be detected using the QUANTOFIX® Total Iron 100 (REF 91344), QUANTOFIX® Total Iron 1000 (REF 91330) and QUANTOFIX® Copper (REF 91304).

Disposal:

Information regarding disposal can be found in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4–30 °C). If correctly stored, the test kit may be used until the use-by-date printed on the packaging.

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