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Rev. 10/2025

Overview

The test is suitable for determining the acute toxicity of freeze-dried luminous bacteria in accordance with DIN EN ISO 11348-3:2023-12.

The test is suitable for surface water, groundwater, seepage water, pore water, sea water, brackish water, wastewater, and aqueous extracts and eluates from sediments.

- Measuring range: 0–100 % inhibition
- Reporting of results:
% inhibition of light intensity in the sample compared to an uninhibited control
 G_L (according to DIN EN ISO 11348 2: LID): Reciprocal of the first dilution level of a sample in which the inhibition of light intensity is less than 20 %
 EC_{20}/EC_{50} : Concentration of a sample that causes an inhibition of light intensity of exactly 20 % / 50 %
- Number of determinations: 100
- Shelf life: 24 months
- Storage temperature:
-21 °C to -19 °C (BioFix® Lumi luminous bacteria)
2 °C to 8 °C (BioFix® Lumi reactivation solution/control solution)

Method

Determination of acute toxicity on freeze-dried luminous bacteria. The measured variable is the natural light output (luminescence) of the microorganism *Allivibrio fischeri* NRRL B-11177 used. The inhibition of light intensity by the sample is determined in comparison to an uninhibited control sample.

Content

- 10 vials BioFix® Lumi Luminous bacteria
- 75 mL BioFix® Lumi Reactivation solution
- 20 mL BioFix® Lumi Control solution

Accessories

- Luminometer (z. B. BioFix® Lumi-10, REF 940008), with suitable cuvettes (e.g. REF 916912) and cuvette holders (e.g. REF 945013)
- Heating block / water bath for temperature control at 15 °C
- Piston pipette 1–5 mL (REF 916909) with pipette tips (REF 916916)
- Piston pipette 100–1000 µL (REF 91677) with pipette tips (REF 91676)
- BioFix® Lumi Osmotic pressure adjustment solution (REF 945602)

Quality control

The enclosed test certificate guarantees that the required validity criteria of DIN EN ISO 11348-3 are met.

The sensitivity of the luminous bacteria must also be checked in your own laboratory according to the standard. The test certificate provides the necessary information, such as standard substances and test concentrations.

The enclosed NaCl-based BioFix® Lumi control solution is used to produce a control sample free of inhibitory substances.

Sampling and preparation

See DIN EN ISO 5667 3 A21.

Turbid sediment samples should be centrifuged or filtered using CHROMAFIL® PES (REF 729242).

Check the oxygen content, which should be ≥ 3 mg/L.

Adjust the pH to 6.0–8.5 using a NaOH or HCl solution.

Check the salt content and adjust to 2 % NaCl. Use one part BioFix® Lumi solution (REF 945062) for adjusting the osmotic pressure to ten parts sample to do this.

If necessary, perform dilution steps (see DIN EN ISO 11348-3, Annex B for details).

Procedure

Reactivation

- Add 6 mL cooled reactivation solution to 1 vial of lumious bacteria
- Mix by shaking
- Stabilize for 5 minutes at 2–8 °C

Dosage

- Add 0.5 mL of luminous bacteria solution to each cuvette
- Incubate for 10 min at 15 °C

Measurement (see the manual for the luminometer used for details)

- Measure initial luminescence I_0
- Add 0.5 mL sample / control
- Mix by swirling
- Incubate for t min at 15 °C
- Measure final luminescence I_t

Interferences

According to DIN EN ISO 11348 3, no pH correction is necessary if the pH value of the sample is between 6.0 and 8.5. pH values < 6.0 or > 8.5 can lead to pH-related inhibition of luminescence.

Undissolved or volatile substances can distort the test result. Turbidity in the sample can cause light loss due to light scattering.

Samples with high oxygen consumption (e.g., samples containing chlorine) or low oxygen content (< 3 mg/L) may show toxicity-independent inhibition of light emission.

A salt concentration of > 35 g/L NaCl or other substances of equal osmolality should not be exceeded, as hyperosmotic effects may occur (see DIN EN ISO 11348 3, Annex D).

Storage

The test is stable until the specified expiration date if stored under the specified conditions. Only thaw the luminous bacteria immediately before reactivation.

Reactivated luminous bacteria must be used within 4 hours.

Reactivated luminous bacteria can be stored at 2–8 °C, although this may result in a decrease in light intensity and sensitivity. Refreezing reactivated luminous bacteria is not recommended.

Disposal

No hazardous substances that are subject to labeling requirements are contained.

Luminous bacteria and test preparations can be disposed of down the drain. However, care must be taken to ensure that toxic samples are disposed of properly.

Information on hazards can be found on the outer label and in the safety data sheet. The safety data sheet can be downloaded at www.mn-net.com/SDS.

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MACHERY-NAGEL GmbH & Co. KG · Valencienner Str. 11 · 52355 Düren · Germany