MACHEREY-NAGEL

QUANTOFIX[®] Relax

- Bedienungsanleitung
- User manual
- Mode d'emploi
- Manual del usuario

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User manual 47

Mode d’emploi 89

Manual del usuario 135

MACHEREY-NAGEL

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QUANTOFIX[®] Relax


User manual

1. Quickstart	51
2. Introduction	53
3. Labels on the device and the box	54
4. Explanation of symbols	55
5. Safety warnings.....	56
6. Technical description	57
6.1 Measuring principle	57
6.2 Functional principle QUANTOFIX® Relax.....	57
6.3 Technical data	58
6.4 CE Certification	58
7. Unpacking and set up	59
7.1 Scope of delivery	59
7.2 Note on ambient surrounding	59
7.3 Setting up the instrument	59
7.4 Description of instrument parts	60
7.5 How to plug the instrument in	61
7.6 How to start the instrument	62
7.7 How to switch off the instrument	62
7.8 How to load the printer paper	62
7.9 How to install batteries (optional)	63
7.10 Instrument self test	63
7.11 Calibration	63
7.12 Use of the instrument	63
7.12.1 Buttons	63
7.12.2 Scroll Menu	64
7.12.3 Round buttons	64
8. Analysis of test strips	65
8.1 Parameter selection	65
8.2 Entering a sample ID	66
8.3 Measurement procedure	66
8.4 Measurement errors	67
8.5 Changing the sequence number ("SN")	67
9. Enter the main menu	68
10. Recall results	69
10.1 Memory List	69
10.2 How to find specific results (filtering)	70
10.2.1 Selecting the date.....	70
10.2.2 Selecting search criteria	70
10.3 Test	71
10.3.1 Display suitable matches.....	71
10.4 How to delete results from memory	72
11. Quality control	73
12. Equipment Settings	74

12.1 Test strips	74
12.2 Favorites	74
12.3 Confidence interval	75
12.4 Correction.....	76
12.5 Password	76
12.6 Printer	76
12.7 Sound	76
12.8 Autostart	77
12.9 Battery	77
12.10 Language	77
12.11 Time/Date	77
12.12 Interface	78
12.12.1 Connecting external devices.....	78
12.12.2 Barcode scanner and keyboard	78
12.13 Customization	78
12.14 Reaction time.....	78
12.15 Print Settings	79
12.16 Data Export	79
13. Maintenance and cleaning.....	80
13.1 How to clean the housing	80
13.2 How to clean the strip holder.....	81
14. Service menu	82
14.1 How to reset the system (Load Default)	82
14.2 How to update the instrument (Program update)	82
14.3 How to run a control strip measurement	83
15. Errors, Sources and Solutions	84
16. Warranty, liability and complaints	85
16.1 General usage comments	85
17. Additional information.....	86
17.1 Electromagnetic Compatibility	86
17.2 Disposal	86
17.3 CAN ICES.....	86
18. Spare parts, Accessories and Consumables	87
19. Manufacturer information	88
19.1 UKCA Authorized Representative	88

1. Quickstart

Unpack the instrument and place it on an even, hard surface. Connect the power supply and turn the equipment on with the On/Off-switch (Pic. 8-⑧). After the self test the start screen will appear on the display.

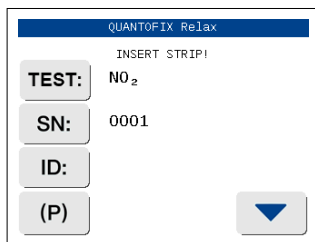Display 1: Start menu

- Pressing **TEST:**, you can choose the parameter to be measured. Once a parameter has been activated, use the test strip as described in the instructions.
- Place the strip on the strip holder (Pic. 1)

Pic. 1: Correct

Pic. 2: Wrong


- Slide or push the strip to the end of the strip holder. Do not touch the reagent pads on the test strip.

The instrument will automatically detect an applied strip. The measurement cycle will be started. A progress bar on the display shows the remaining analysis time.

NOTICE

If "Autostart" (Chapter 12.8 "Autostart") is deactivated, the measurement must be started using the start control panel .

At the end of the measurement the result will be displayed on the screen and transferred to the printer and interfaces.

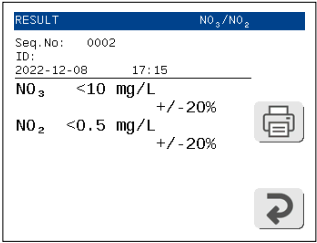Display 2: Result

By pressing the printer symbol  the result can be printed again. Choosing the return panel  will lead back to the start screen.

Another analysis may be started by applying the next test strip.

NOTICE

To start a new measurement it is not necessary to go back to the start screen. A new strip is detected at any time and the measurement is then started automatically.

2. Introduction

MACHEREY-NAGEL

The QUANTOFIX® Relax is a reflectometer for the evaluation of test strips from MACHEREY-NAGEL. Only use test strips from MACHEREY-NAGEL that are suitable for instrumental evaluation.

The measurements are carried out under standardized conditions. Results may be displayed, printed and sent to a computer. Instrumental evaluation eliminates potential sources of error inherent to visual evaluation. Results will not vary due to subjective color perception or different light sources. Additionally, errors resulting from manual result administration are avoided. Featuring a clear touch screen and intuitive user guidance, the QUANTOFIX® Relax is ideal for quick, efficient and comfortable work.

3. Labels on the device and the box

Pic. 3: Exemplary carton label

Pic. 4: Exemplary type plate label

4. Explanation of symbols

The device and its packaging contain the symbols and terms listed below with the following meanings:

This symbol indicates that there is a risk of electric shock.

This symbol indicates that the designated area can become hot and should not be touched without suitable protective measures.

This symbol indicates that the chemicals used have a corrosive effect. Observe the safety measures in the laboratory and wear the prescribed protective equipment. Observe the instructions in the current safety data sheets (SDS) of the products used.

This symbol indicates that there may be a danger by using flammable substances.

Caution

USB interface

I/O

On / Off

LAN-interface

RS232

COM-Interface

Direct current

Keep dry

Temperature limits

Consult instructions for use or consult electronic instructions for use

Do not dispose of the device in common household waste

Manufacturer

Catalogue number

Serial number

Date of Manufacture

CE Mark

UKCA Mark

5. Safety warnings

BE SURE TO READ THE FOLLOWING SAFETY INSTRUCTIONS CAREFULLY BEFORE USING THE DEVICE.

Failure to follow these instructions may result in serious injury to the operator, malfunction or damage of the device.

Keep this manual in a safe place for future reference.

Follow the safety notes and instructions in the operating manual and observe the stickers and notices on the device.

In case of a test strip sled blockage, switch off the device and turn it back on after 10 seconds. If the blockage persists, do not continue using the device and contact technical support.

If liquids have entered the interior of the device, do not use it any further and contact technical support.

If there are any cracks or holes in the housing, stop using the device and contact technical support.

If the software freezes, switch off the device and turn it back on after 10 seconds. If the software remains unresponsive, do not continue using the device and contact technical support.

Do not work on internal parts of the unit. Non-compliance will invalidate any warranty claim.

The following safety warnings are used throughout the manual:

DANGER

Indicates an imminent or potentially hazardous situation which, if not avoided, will result in death or serious injury.

WARNING

Indicates an imminent or potentially hazardous situation which, if not avoided, could result in death or serious injury.

CAUTION

Indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury.


NOTICE

Indicates a situation which, if not avoided, may result in malfunction or damage of the instrument. Information that clarifies facts in the text and that requires special attention.

6. Technical description

6.1 Measuring principle

The test strip moves below a fixed measuring head on a sled with an embedded reference pad. The reflectometric analysis of the test strip and the reference field take place during withdrawal and release of the sled.

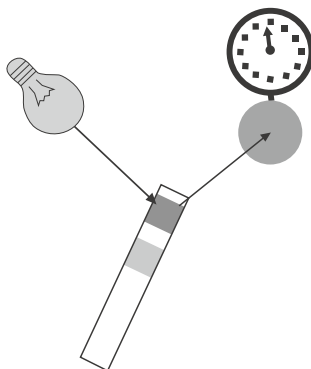Pic. 5: Measuring Principle

The strip is illuminated with an LED and a detector registers the intensity of light reflected by the test strip at three different wavelengths. Using an internal calibration, the intensity of the reflected light is correlated to the concentration of the analyte.

6.2 Functional principle QUANTOFIX® Relax

A measurement is started by placing a strip on the holder. If the Autostart-feature is turned off, the measurement is started by pressing the start panel  in the display. The result is shown on the display, printed out and released via the interfaces after the measurement has been completed. After three minutes the instrument will go to stand-by. Touching the screen will reactivate the instrument.

All user inputs are performed via the touch-screen (Chapter 7.12 “Use of the instrument”).

6.3 Technical data

Specification	Description
Size	7.5 cm × 16 cm × 20 cm (H × W × D)
Weight	710 g (without batteries and power pack)
Power supply	External power pack 100–240 V, Output 9 V, 1.5 A Alternative: battery operation with 6 mignon batteries 1.5 V (AA)
Operating conditions	10 °C–40 °C; 20 %–80 % rel. humidity (non-condensing), up to 3000 m
Pollution degree	2
Protection Class	Device: III Power supply: I Similar to UL 50E*: Type 1*
Overvoltage category	Device: II Power supply: I
Transport condition	-10 °C–45 °C, max. 80 % humidity (non-condensing)
Storage condition	10 °C–40 °C, max. 80 % humidity (non-condensing)
Application Area	Commercial indoor laboratory

* This classification for the enclosure corresponds to typical Type 1 enclosures in terms of design and protective effect, but it has not been tested in accordance with UL 50/UL 50E and is not UL-certified.

The maximum sound level and maximum sound pressure are subject to the limits specified in ISO 3746 or ISO 9614-1.

6.4 CE Certification

The CE mark declares that the product complies with the harmonization legislation of the European Community listed below:

European Directive 2011/65/EU on the restriction of the use of certain hazardous substances in electrical and electronic equipment (RoHS 2)

This device complies with the following regulations:

IEC 61010-1

IEC 61326-1

FCC47 CFR Part 15B, ICES-003 Issue 6

ISED ICES-001 Issue 5 - CAN/CSA CISPR 11:19

7. Unpacking and set up

7.1 Scope of delivery

- ① QUANTOFIX® Relax reflectometer
- ② Power pack 100 – 240 V, 47/63 Hz, 9 V, incl. 4 adapters
- ③ Printer paper
- ④ User manual (this booklet)
- ⑤ Quick start guide (different languages)

Pic. 6: Content

Check all delivered parts for visible damages. In case a part is damaged, please contact your local distributor or MACHEREY-NAGEL.

Read the operating manual for QUANTOFIX® Relax carefully before the first startup in order to ensure an error free operation.

NOTICE

Position the device with at least 5 cm of free space at the back to ensure safe access to the power switch.

7.2 Note on ambient surrounding

If the device is exposed to higher temperature fluctuations (e.g. after transport or distribution), it must be switched on not before sufficient acclimatization is given. The device should not be used close to electrical fields (e.g. by microwaves, radio units et cetera). In worst case the measurement results can be affected.

7.3 Setting up the instrument

Place the instrument on a hard, even surface where humidity and temperature are fairly constant. Make sure that the instrument is allowed to acclimate to room temperature prior to use.

Make sure that you

- Do not place the instrument near strong electromagnetic fields
- Do not place the instrument near heating plates, ovens or radiators
- Do not expose the instrument to strong light sources (i.e. direct sunlight)

7.4 Description of instrument parts

CAUTION

Do only connect the intended devices to the corresponding interface. If you connect another device to the corresponding interface, the analyzer or the connected device may be damaged. Check all cables prior to use and verify the proper connection.

CAUTION

Multiple hazards: Only qualified personnel should carry out instructions described in this step.

WARNING

Risk of electric shock: Make sure that the power cord is not damaged. Check the suitability of the power source for the equipment.

CAUTION

Risk of injury due to the danger of tipping over due to slipping. Place the device on a flat surface. Do not stack the device.

WARNING

Fire hazard: Take care not to overload the mains socket. There is a risk of overload and fire.

Make sure that the power cord is not damaged. Check that the power source used is suitable for the equipment.

WARNING

Risk of injury: Familiarize yourself with the device before working with it and read this document carefully. Do not use the device unless you have received instruction in its use.

NOTICE

Defects of the power supply and housing can lead to malfunction of the device. If the device shows an apparent breakage of the housing or a damaged power supply, it must be taken out of operation.

Defects of the power supply and housing can lead to malfunction of the device. If the device shows an apparent breakage of the housing or a damaged power supply, it must be taken out of operation.

Pic. 7: Front view

Pic. 8: Backside view

Actuator	Function
1. Touch-Screen	Control of equipment functions
2. Test Strip Slide	Test strip retainer and autonomous start of analysis
3. Printer Flap	Opening the printer flap for paper replacement
4. USB A -Interface	Connection of a keyboard or a bar code scanner ²
5. USB B - Interface	Connection of a computer (cable length ¹ max. 3 m)
6. LAN - Interface	Currently not supported
7. Serial Interface	Connection of a computer (cable length ¹ max. 3 m)
8. On/Off Switch (I/O)	Turning the equipment on and off
9. Mains Connection	Contact for the provided power pack

¹ We recommend to use shielded cables.

² We recommend to make use of the hand scanner from MACHEREY-NAGEL

7.5 How to plug the instrument in

NOTICE

Only the associated power supply (REF 930995) may be used.

Pic. 9: Power pack

Pic. 10: DC in

Four adapters are provided for adapting the power pack to the available mains connection. The connector pins sealing is removed and the adapter matching the mains connection is

plugged on to the power pack (Pic. 9). After plugging the power pack cable into the jack "DC IN" (Pic. 8-g) and connecting the power pack to the power socket the QUANTOFIX® Relax is ready for operation.

To ensure safe functioning of the instrument, only use the included power supply!

7.6 How to start the instrument

You can find the On/Off switch on the back of the instrument. Turn on the instrument by pressing the switch. Right after switching on, the instrument performs a self-test. The touch screen lights up and the start screen appears. If a serious error is detected during the self-test, no measurements can be performed. In this case, contact the MACHEREY-NAGEL technical service.

7.7 How to switch off the instrument

Always switch off the device after use by pressing the on/off switch. If you do not use the device for a longer period of time, disconnect the device from the power outlet.

7.8 How to load the printer paper

Pic. 11: Printer A

Pic. 12: Printer B

Open the printer flap by pressing the rectangular key next to the printer flap (Pic. 11).

Pic. 13: Printer C

Pic. 14: Printer D

Unroll the paper roll by 5 cm and place the roll in the paper compartment with the end on the lower side. Fix the end of the paper to the housing with your finger while closing the flap (Pic. 13 + Pic. 14).

7.9 How to install batteries (optional)

The QUANTOFIX® Relax can be operated with 6 type AA batteries independent of the mains supply. The battery compartment is on the underside of the device. Notice the designated polarity (+/-) marked on the battery compartment while inserting the batteries.

Pic. 15: Battery compartment

7.10 Instrument self test

The instrument will perform an automatic self test each time it is turned on. If an error message appears, the instrument will not start measurements. In this case, please contact your local service provider.

7.11 Calibration

The instrument will perform an automatic calibration each time a test is performed.

7.12 Use of the instrument

All user inputs are done via a touch-screen (touch-display). All functions are activated directly by slight pressure with the finger on explicit pictograms or text representing the menu items.

7.12.1 Buttons


Framed areas react to pressure and trigger the action linked to it. The caption of an area describes its function.

Examples:

- | | |
|--|--------------------------|
|  | Confirm / perform action |
|  | Cancel action |
|  | Check mode |
|  | Memory |
|  | Enter main menu |

7.12.2 Scroll Menu

Press the up-and-down arrows on the right side of the screen to scroll through a list of information on the left side of the screen. The desired information on the left side is highlighted.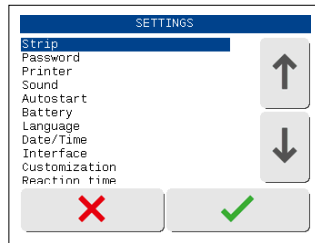


Display 3: Selective lists

Pressing  will select the highlighted line. You can leave the menu by pressing .

7.12.3 Round buttons

These buttons typically appear on screens that require a selection among serial items. The button with a filled circle is the current selection.

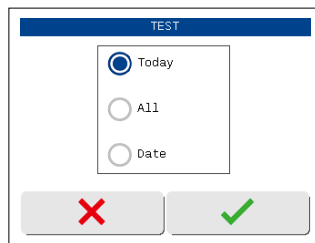Pressing the circle will activate a selection. Save your selection by pressing . Pressing  will quit the menu without performing any changes. Pressing the entry of the filled circle will deactivate the function in some menus Chapter 12.12 "Interface"

8. Analysis of test strips

The instrument is designed for the quick and easy measurement of test strips.

The following points outline set-up options with regards to measurements as well as the actual measurement procedure.

WARNING

Possible dangers of contact with chemical and biological substances.

Working with samples, reagents and corresponding accessories is associated with dangers.

Wear suitable protective equipment when working with test kits. Observe the safety data sheet (SDS) of the test kit used.

8.1 Parameter selection

NOTICE

Make sure to remove excess liquid by blotting the test strip carefully on a lint-free cloth.

NOTICE

If the autostart mode (Chapter 12.8 “Autostart”) is deactivated, the analysis needs to be started by pressing  on the touch-screen.

The instrument cannot detect different test strips for different parameters or ranges automatically. Therefore, initially you have to activate the test strip you want to measure. The currently activated test is displayed in the start screen next to the button **TEST:**. Always make sure that the test strip you are using corresponds to the activated test.

To activate a particular test for a measurement, press **TEST:**. A new screen appears which includes up to 5 pre-defined tests (see Chapter 12.2 “Favorites”).

To activate one of these favorites, press the corresponding button. The start screen appears and the desired test is activated and is displayed next to the **TEST:** button.

To activate a test that is not part of the favorites list, press . A list with all available tests appears. Use the arrows to select a test. Press  to confirm your selection, the start screen appears and the desired test is activated and displayed next to the **TEST:** button. If you leave the list by pressing , your selection is not adopted.


The instrument saves the last used test and keeps it activated in the start screen.

After the measurement, the instrument will release the analyzed test strip which can now be discarded. The result is displayed on the screen and is transferred via the interfaces and/or printed according to equipment settings.


For additional information on the test strip, please read the instruction leaflet that comes with the strips.

8.2 Entering a sample ID

ID: shows the sample ID for the next measurement. To enter or change the sample ID, press **ID:**. The input menu appears and you can enter the desired sample ID. The following measurements are then always run and saved using the same sample ID. Using the search mode you can retrieve the results from the instrument memory. You can also enter the sample ID using a barcode scanner or a keyboard.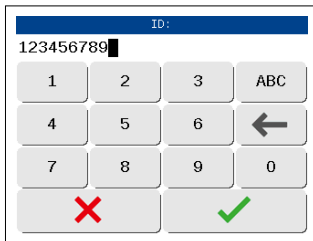

Display 4: Entering ID (numeric)

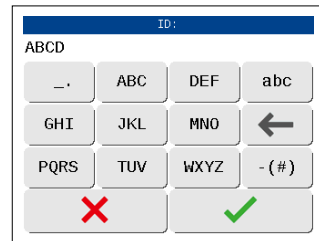Display 5: Entering ID (letter)

- Using a standard PC-keyboard: Connect the keyboard to the USB-A interface in the backside of the instrument.
- Using a bar code reader: Connect the barcode reader to the USB-A interface in the backside of the instrument.

8.3 Measurement procedure

NOTICE

It is important to place the test strip at the far end of the slide's recess. You can easily move the strip forward on the recess using your finger (make sure to not touch the test field!).


NOTICE

The reaction time is dependant on the test and programmed into the instrument. Therefore, different tests have different waiting times. In general though, the time to wait is hardly ever longer than 60 seconds.

If the autostart function is disabled, you have to start a measurement pressing  (Chapter 12.8 "Autostart").

Measuring test strips with the QUANTOFIX® Relax is very easy. To start a measurement, simply place the test strip on the test strip slide with the test pad facing upwards and forward. The end of the test strip has to touch the end of the slide's recess. A sensor (little gap within the slide) detects the test strips and starts the measurement automatically.

A progress bar appears showing the remaining reaction time. After a few seconds, the test strip is drawn into the instrument's housing. After the measurement, the slide moves out of the housing again and the result is displayed in the screen. The measured test strip can be discarded into the domestic waste.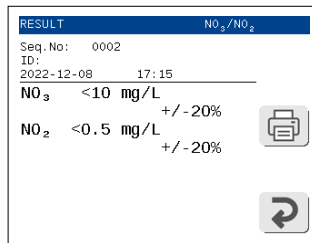

Display 6: Result

8.4 Measurement errors


If the display shows a measurement error (e.g wrong strip) instead of a result please read the instructions in Chapter 15 "Errors, Sources and Solutions".

Repeat the measurement. In case of permanent errors please contact MACHEREY-NAGEL service.

Measurement errors will be printed, send and stored according to the instrument settings.

8.5 Changing the sequence number ("SN")

Pressing  in the start menu brings up a numerical pad. Enter a new sequence number using the keys on the pad. All following measurements will now be counted starting from this number. For every new day, the sequence number is set back to 0001.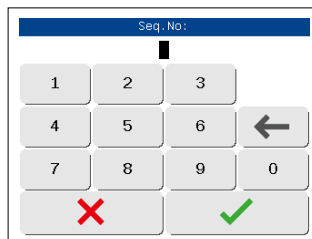


Display 7: Seq.-Input


9. Enter the main menu

Pressing  on the start screen will bring up the main menu. Within the main menu there are three options.

-  allows you to display, search and delete stored data.
-  enables quality control measurements with special MN control strips.
-  gets you to the instrument settings.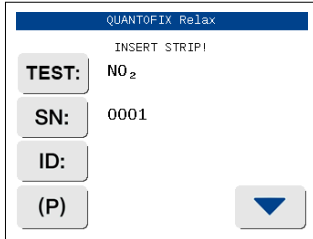

Display 8: Start menu

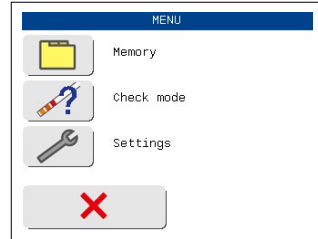
Display 9: Main menu


10. Recall results

The QUANTOFIX® Relax has an internal memory for storage and recall of measurement results. Every result is automatically saved after the analysis. When the memory is full, new data will overwrite the eldest saved dataset.

Access the memory by pressing  in the main menu .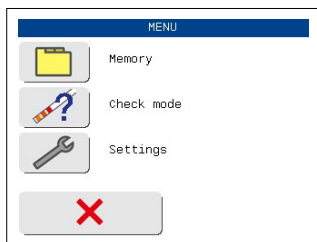

Display 10: Main menu

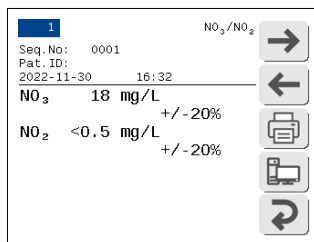
Display 11: Memory mode

Within the memory menu, you have three different options. You can display a complete list with all data sets, search for specific data sets or delete the entire memory.

10.1 Memory List

Pressing  to display all data sets in the memory. A new screen appears, showing the result for the latest measurement. The results are always sorted according to the measurement date. The heading shows the number of the currently displayed data sets and the total number of data sets within the memory.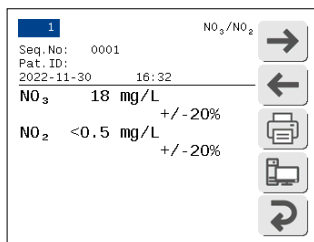

Display 12: Memory contents


For the respective result, sequence number, sample ID, measurement date/time, test and the actual result are displayed. On the bottom side of the screen are 5 activity buttons.

Press  and  to choose between different data sets. Additionally, you may press  to print the displayed data set or press  to send the data to a PC.

Press  to get to the memory menu.

10.2 How to find specific results (filtering)

Press  to search the memory for specific results. To do so, you can freely combine three search criteria (test, date, parameter). To start a search, all criteria need to be defined.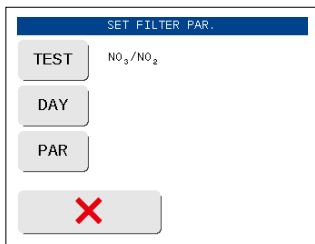


Display 13 shows the 'SET FILTER PAR.' menu. It has a title bar 'SET FILTER PAR.' and three buttons: 'TEST' (with 'NO₃/NO₂' next to it), 'DAY', and 'PAR'. At the bottom is a large button with a red 'X'.

Display 13: Filtering

10.2.1 Selecting the date

By pressing  you will reach the menu displayed below (Display 14).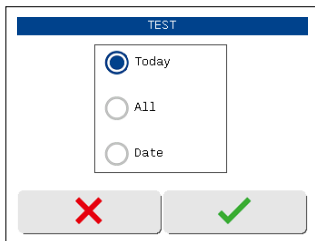


Display 14 shows the 'TEST' menu. It has a title bar 'TEST' and three radio buttons: 'Today' (selected), 'All', and 'Date'. At the bottom are two buttons: one with a red 'X' and one with a green checkmark.

Display 14: Select day

Set the day with the buttons. Selecting "Date" will bring up a screen with the list of available dates (only days with measurements are shown on the screen). Select the desired date with the up-and-down arrows and confirm your selection by pressing . After confirmation your selection will be displayed on the screen "SET FILTER PAR."

10.2.2 Selecting search criteria

Pressing  in Display 13 brings up Display 15.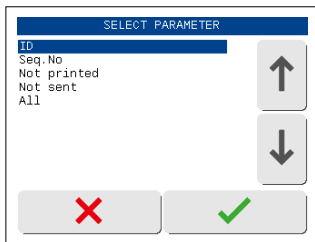


Display 15 shows the 'SELECT PARAMETER' menu. It has a title bar 'SELECT PARAMETER' and a list of parameters: 'ID', 'Seq.No', 'Not printed', 'Not sent', and 'All'. To the right of the list are two arrow buttons (up and down). At the bottom are two buttons: one with a red 'X' and one with a green checkmark.

Display 15: Select parameter

Use the arrow keys to select the desired criteria and confirm with . The filter criteria will be displayed on the filter settings screen (Display 17).

10.3 Test

Press **TEST** to determine the type of test strip that generated the result you are looking for. Once you press the activity button, a new screen with a list of all available tests appears.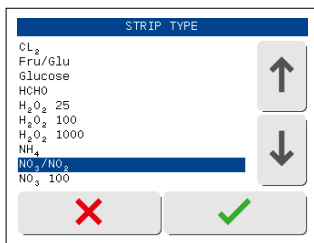


Display 16: Test selection

Use the arrow buttons to choose the test you are looking for. Press  to confirm your selection. Your selection will appear in the search menu. If you leave the list by pressing , your selection is not adopted.


10.3.1 Display suitable matches

After setting "Day" and "Parameter" the search may be started by pressing  (Display 17).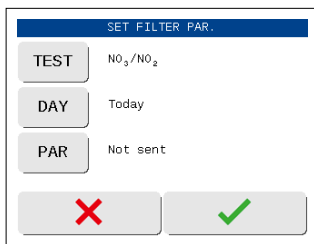

Display 17: Filtering

When suitable matches are found, an option will appear, allowing to print the datasets , send them to a PC , or display them on the screen  (Display 18).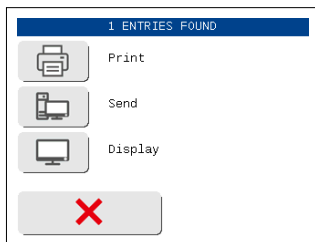

Display 18: Search result

If no matching results are found, the equipment returns to the memory menu.

10.4 How to delete results from memory

Pressing  will delete all data in the memory. You need to confirm this again on a further screen. The quality control measurements are not affected by this action.


11. Quality control

The instrument performs a self-calibration whenever it is turned on. Therefore, in general, the instrument does not need to be calibrated. To check instrument performance, we recommend to run multiple standard measurements with known concentrations within the tests range. If such results are not within the test's confidence interval, (see Chapter 12.3 "Confidence interval") please contact your distributor or MACHEREY-NAGEL directly (see Chapter 19 "Manufacturer information").

For additional quality control assessments, you can also use the instrument's check mode and special control strips (see Chapter 14.3 "How to run a control strip measurement").

12. Equipment Settings

Enter the main menu by pressing the menu key  to reach the "SETTINGS" display press ). Use the arrow buttons to choose an entry and confirm your choice with .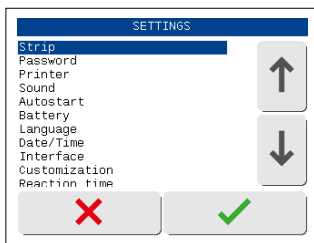

Display 19: Admin settings

12.1 Test strips

Within the "Test strip" menu there are three more options. Favorites allow you to define test strips for the "Favorites" menu (see Chapter 12.2 "Favorites"). Additionally, you can activate or disable the confidence interval message in the display.

The option "Correction" allows for the adjustment of the factory-provided parameter calibration to account for matrix effects and interferences during the analysis. Use the arrow buttons to choose an entry and confirm your choice with .


If you press  you get back to the settings.

12.2 Favorites

You can define up to 5 different test strips as favorites. These favorites are saved and will appear in the "Favorites" menu. You can access the favorite's selection menu via **TEST:** in the start screen. The favorites option allows you to pre-define your most frequently used tests, so that you can quickly and easily switch between test strips. The menu shows a list with all tests and parameters that are programmed for evaluation in the instrument. On the left side of the

screen, you can see the test names. On the right hand side of the name, directly next to the arrow buttons, you can see a confirmation symbol, given the respective test has been marked as a favorite in the first place.

To add a test to the favorites list, use the arrow keys to navigate to the desired test and then press the Favorites icon (star) located between the arrow keys. Confirm your selection by pressing the confirmation button .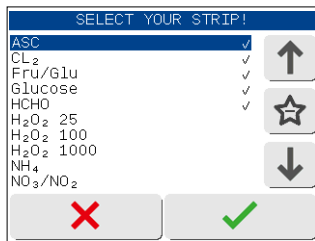

Display 20: Menu for selecting favorites

The confirmation symbol appears behind the test, it has been included into the favorite menu. You can deactivate parameters from the favorite list in the same way.

The favorite list can include only 5 tests. If there are already 5 tests activated within the list, you have to deactivate a test first in order to choose a new favorite.

To leave the menu, press , and you will get back to the start screen.

12.3 Confidence interval

Naturally, test strip results are subject to slight deviations even when they are generated photometrically. For every test there is a percentage based confidence interval in which the true value can be found. The instrument can optionally display the confidence interval (in +/- percent) underneath the result after every measurement. To activate or deactivate the confidence interval, respectively, choose the desired choice button from the option field and confirm with .

If you press , none of your selections is adopted.

12.4 Correction

WARNING

Faulty measurement results due to manipulated calibration are the sole responsibility of the operator of the equipment!

The correction mode allows for the adjustment of the MACHEREY-NAGEL parameter calibration and affects the result in terms of an additional calculation step. The menu offers the opportunity of correcting the result with a linear correction function.

By selecting the correction menu four values appear, that can be edited by choosing each of them. Selecting this function always affects the parameter displayed in the strip menu (see Chapter 12.1 “Test strips”).

The meaning of the values A to D is as follows:

A = gradient of the correction function

B = y-axis intercept of the correction function

C = new lower measurement limit

D = new upper measurement limit

Confirming your entry with  stores the settings. If you press , none of your selections is adopted. By pressing the “Reset”-button the values will be set back to factory settings.

Once you select one of these values, the input menu appears and you can enter a value (Restrictions: The value for A must be > 0 . Value D must be bigger than value C, if values C and D are equal to zero, the factory-provided measuring range will be adjusted depending upon the entered values for A and B).

Confirming your entry with  stores the value. If you press , the former value will be kept.

The modification of a parameter will be indicated with a star in the parameter list display as well as on the result display and printout. Also the result in the memory will contain this indication, which cannot be removed by resetting the values to factory-settings.

12.5 Password

You can protect access to the instrument settings using a password, thus restricting the number of user who can access the settings. To install a password, press “On” in the option field and confirm with . If you press , none of your selections is adopted.

The input menu appears and you can enter a password. Once you leave the settings and then go back again, the instrument requires the respective password.

12.6 Printer

In the printer menu, you can disable or enable the printer. Press the respective choice button in the option field and confirm with . If you press , none of your selections is adopted.

12.7 Sound

In the “sound” menu, you can disable or enable the instrument sound. Press the respective choice button in the option field and confirm with . If you press , none of your selections is adopted.

12.8 Autostart

The instrument has an autostart function which detects a test strip on the slide and automatically starts the measurement process.

In the autostart menu, you can disable or enable the autostart. Press the respective choice button in the option field and confirm with . If you press , none of your selections is adopted.

If the autostart function is disabled, the start screen shows . In this case you have to press the activity button to start a measurement.

12.9 Battery

You can make some adjustments to conserve energy. Within the menu you will find to option fields. Here you can enable or disable the screen's backlight function and the printer. Press the respective choice button(s) and confirm with . If you press , none of your selections is adopted.

12.10 Language

The instrument can be set to 15 different languages. In the "Language" menu, you can choose between the different languages. The menu contains a list of all languages available. To activate a particular language, select the language using the arrow buttons and confirm with . The instrument changes the language automatically and goes back to the settings menu. If you press , none of your selections is adopted.

12.11 Time/Date

Select "Date/Time" in the setting menu. To change, press on the respective number. A numerical pad appears. Enter the correct number and confirm by pressing .


The date may be formatted in three ways. The active format is shown on a button  on the right hand side of the date (Display 21). Select a date format by pressing this button.

Displayed Format Abbreviation	Meaning	Example
YMD	Year - Month - Day	2007-12-17
DMY	Day . Month . Year	17.12.2007
MDY	Month / Day / Year	12/17/2007

Changing the displayed format will influence the time format during the use of the instrument and the printout, but will not affect the format of the data send via the interfaces.

The time format may be changed to 12 or 24 hours with the button displayed next to the time

24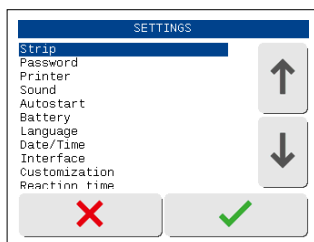

Display 21: Date / time

12.12 Interface

In the "Interface" menu, you can disable or enable the instrument's interface.

The instrument can send data only if the interface is enabled. Press the respective choice button in the option field and confirm with . If you press , none of your selections is adopted.

12.12.1 Connecting external devices

The instrument can be connected to a barcode scanner and/or keyboard as well as to a computer. If connected to a computer, the instrument can only send, but not receive data (unidirectional communication).

12.12.2 Barcode scanner and keyboard

Both barcode scanner and keyboard can be connected via USB-A. Using a barcode scanner (Chapter "18. Spare parts, Accessories and Consumables") or keyboard enables you to enter sample IDs or to make entries in open input fields.

12.12.3 Transferring data to a PC

You can connect the instrument to a computer using its USB or RS/232 interface (Chapter 12.12 "Interface"). For the transfer of stored data from the device to a computer, the data export software for strip reader from MACHEREY-NAGEL is available free of charge. The data received by the data export software for strip reader can be stored either as a Microsoft® Excel / Open Office table or as a text file (.txt).

For more information about the data export software, please refer to the following website: www.mn-net.com/quantofix-relax-geraete-update

For further information regarding connection to computers or laboratory information systems, please contact your distributor or MACHEREY-NAGEL directly (Chapter 19 "Manufacturer information").

12.13 Customization

Select "Customization" in the settings menu. The first two lines of the printout may be filled with a user-specific identifier. Each line contains 23 characters.

To enter the text an external keyboard or the alphanumerical pad on the display can be used. The keys on the touch-screen are linked to several letters. Repeated pressing within half a second switches through the letters displayed on the key.

12.14 Reaction time

In the "Reaction time" menu, you can decide whether the instrument will wait for the strips reaction time before starting a measurement. By default, the reaction time is enabled. The reaction time is then shown as a progress bar once the strip is placed on the strip holder.

If you disable the reaction time, the instrument draws the strip into the housing immediately and starts the measurement irrespective of the strip's individual reaction time. Hence, you need to place the strip onto the strip holder as soon as the reaction time has run out. In such a way, you can potentially generate a higher strip throughput.

To activate / deactivate the reaction time, press the respective choice button in the option field and confirm with . If you press , none of your selections is adopted.


The switched off reaction time will be indicated with a hash on the result display and printout.

12.15 Print Settings

The last menu item print settings allows you to print an overview over the adjustments made settings. Select the entry using the arrow buttons and press . The overview is printed automatically.


12.16 Data Export

To enable data export via a USB flash drive, connect a USB flash drive to the USB A port on the rear of the Relax device and enable data export in the settings.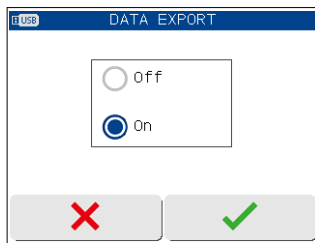

Display 22: Data export settings



Whether a USB flash drive is connected and recognized is indicated by a USB symbol in the upper left corner of the screen.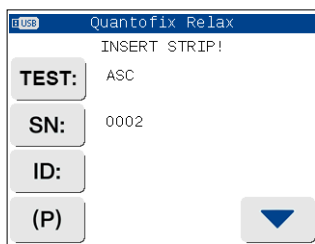

Display 23: Home screen with USB symbol

Measurement data is automatically transferred to the USB flash drive after each completed measurement. The device creates a separate export file for each calendar day on the USB flash drive and continuously adds all measurements recorded on that day to the file.

Alternatively, data can be exported to the USB flash drive at a later time. For this purpose, either individual measurements or the complete device memory can be exported using the USB symbol.

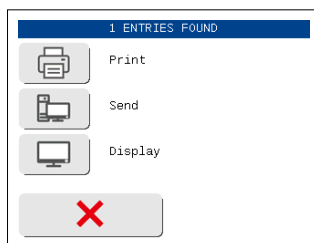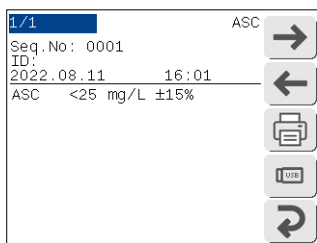Display 24: Single measurement and memory selection in data export mode

13. Maintenance and cleaning

NOTICE

For all cleaning work, the unit must be switched off and disconnected from the power supply.

WARNING

Fire hazard. Do not use flammable detergents or organic solvents to clean the unit.

NOTICE

Do not clean the device with acetone or similar products.

NOTICE

Malfunction or damage of the device due to incorrect maintenance. Maintenance of the device may only be carried out by qualified personnel

NOTICE

Only use accessories and spare parts approved by the manufacturer. If parts not approved by the manufacturer are used, the operator is responsible for the conformity of the device and the warranty will be voided.

To ensure correct functioning of the instrument for a long time, the instrument needs to be cleaned and maintained carefully.

13.1 How to clean the housing

NOTICE

Use commercially available disinfection wipes to clean and disinfect the instrument from the outside.

NOTICE

Do not use liquids that can enter the instrument as these may cause permanent damage of the instrument.

Wipe the outside of the device with a damp cloth, if necessary with a mild cleaner, and then dry the device with a soft cloth. Make sure that no moisture enters the device or the printer. Clean the touchscreen with a lint-free cloth. Avoid scratching the display. Do not use ballpoint pens or other pointed objects to operate the touch screen.

13.2 How to clean the strip holder

Wipe off sample residues from the strip holder with a lint-free cloth after each measurement. This prevents carry-over and drying of sample residues.

To remove the strip holder for more comprehensive cleaning, make sure to turn the instrument off. Pull the strip holder from the metal sled below (Pic. 16). Wipe down the strip holder with disinfection wipes. When removed from the instrument, you may also use water with a soft brush and/or liquid disinfectants.

After cleaning, insert the holder on the metal sled (Pic. 17). The rectangular notches of the transport mechanism and the strip holder must be placed on top of each other (Pic. 18). Use reasonable force to push the strip holder completely onto the metal sled. Doing so the metal sled and the strip holder will be pushed back completely into the housing. If you get an error message after turning the instrument on: Turn the instrument off and push the strip holder again on the metal sled with more force.

Pic. 16: test sled
(bottom view)

Pic. 17: Reinsertion of the
test strip holder

Pic. 18: Check alignment of
notches


NOTICE

Ensure that the strip holder is completely clean and dry before inserting.

14. Service menu

The QUANTOFIX® Relax has a password protected service menu. To enter the service menu press the touch-screen three times during the self test after turning the equipment on (Display 25). Upon request input the PIN "1234".

A selective list with different menu items appears.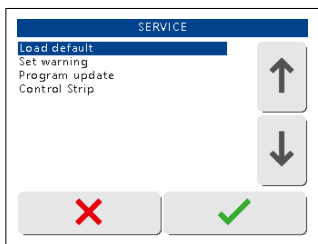

Display 25: Service menu

14.1 How to reset the system (Load Default)

Select "Load default" from the service menu. Confirming this entry will call up a safety question. Pressing  will reset the delivery status.

NOTICE

All settings modified by the user will be cleared!

14.2 How to update the instrument (Program update)


Select "Program update" from the service menu. The instrument now expects the upload of a new firmware. Follow the instructions that come with the update-file to finalize the update process.

14.3 How to run a control strip measurement

NOTICE

Do not touch the test fields of the control strip with your fingers.

The check mode for control strips allows you to check the correct functioning of the instrument using special color control strips. You can order these control strips from MACHEREY-NAGEL. Choose the entry "Control strip" in the service menu. The instrument asks for insertion of the colored control strip.

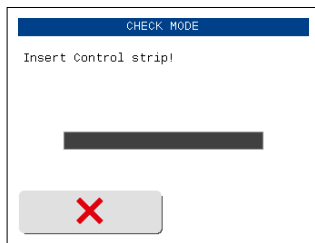Display 26: Control strip

The measurement of the control strip starts automatically as soon as the strip is placed on the sled, if the autostart is activated. Once the measurement has been completed, the measured values are immediately displayed and printed. The result of the measurement is not stored. The values are so called remission values for the pre-colored test pads on the control strip. These values have to be compared to a set of standard values from the control strip package instructions.

For additional information on the control strips or in case the values do not fit the expected value ranges in the control sheet, please consult the package insert, or contact MACHEREY-NAGEL directly.

15. Errors, Sources and Solutions

The instrument differentiates between different types of error messages. Advisory error messages are of less importance and are displayed on the screen during handling the instrument (e.g. Please remove strip). When the corrective action has been performed, the analyzer removes the message from the screen.

If an error message occurs during testing it will be displayed on the screen or in the result view instead of the result. Messages are displayed in plaintext and are self-explanatory.

Error	Source	Solution
"Wrong Position"	The strip hasn't been pushed into the strip retainer far enough	New measurement, place strip in right position
"Printer out of paper"	Paper roll empty or printer flap open	Replace paper and close printer flap
"Battery Low"	Batteries are low	Exchange batteries or use power pack
Instrument doesn't start	Power supply not installed or defect	Check whether all connections are plugged in and whether the power socket is functioning
"Mech. error"	Hardware error	Please check the correct positioning of the strip holder. Please contact your local distributor of MACHEREY-NAGEL directly.
"Measurement error"	An error during the measurement occurred.	Please repeat the measurement.

In case an error cannot be solved by the aid of the instructions above, please contact your local distributor or the MACHEREY-NAGEL Service.

16. Warranty, liability and complaints

The warranty for this equipment has a duration of 24 months from the date of purchase. The original copy of the bill serves as a certificate and must be submitted in case of assertion of a warranty claim. The warranty expires in case of improper handling and/or maintenance of the equipment; it does not comprise defects due to the external power supply.

The warranty is limited to the repair of faulty parts or – at our sole discretion – to the delivery of a faultless substitute. The warranty period of 24 months is not affected by claiming on the warranty during this period. There is no right of withdrawal.

Further claims are excluded. Hereunto we count in particular all claims for damages evolving from consequential damages or indirect damages.

Additionally the relevant version of our general sales terms and delivery conditions apply as printed on all price lists.

16.1 General usage comments

NOTICE

Due to the reasons stated above, MACHEREY-NAGEL declines any responsibility for the correct functioning of the instrument if it has been opened or used improperly and objects to any warranty claims made in such a case. MN's warranty and liability obligations as to product(s) or to any party on account thereof shall be void and of no effect if the product(s) or any part thereof have been subject to unauthorized change, abuse, misuse, accident or alteration, unauthorized application or installation not in compliance with the product(s) specifications as set forth in the product leaflets / instructions hereof ("unauthorized action") and such unauthorized action resulted in the non-conformance or defect of the product(s).

17. Additional information

17.1 Electromagnetic Compatibility

Target environment of device use: Application in the commercial indoor laboratory environment.

- This device has been developed and tested in accordance with CISPR 11 Class A. It can cause radio interference in a domestic environment. In this case, you may need to take action to correct the interference.
- Use of this equipment in a dry environment, especially if synthetic materials (synthetic fiber clothing, carpets, etc.) are present, can cause destructive static discharges that can cause erroneous results.
- Do not use this device near sources of strong electromagnetic radiation, as they may interfere with proper operation.
- This device is in accordance with the applicable parts of the requirements for emission and immunity described in the IEC 61326 series of standards.
- Before use, the electromagnetic environment must be assessed.

EMC Class A according CISPR 11:

The device can be used in all institutions excluding residential areas and areas which are directly connected to the public power supply, if the public power supply supplies buildings for residential purpose.

17.2 Disposal

NOTICE

Disposal using public waste disposal facilities is not permitted. In case of disposal, please contact your MACHEREY-NAGEL representative.

Disposal according to EU Directive 2012/19/EU. In accordance with the EU Directive 2012/19/EU MACHEREY-NAGEL takes back the old device and disposes it free of charge.

! DANGER

Dispose hazardous, infectious or biologically contaminated materials in a safe and acceptable manner and in accordance with all local and regulatory requirements.

17.3 CAN ICES

This equipment is classified as industrial, scientific, and medical (ISM) equipment and complies with the requirements of the Canadian Interference-Causing Equipment Standard ICES-001, issue 5, Class A.

18. Spare parts, Accessories and Consumables

NOTICE

Use only accessories and spare parts approved by the manufacturer. If parts not approved by the manufacturer are used, the operator is responsible for the conformity of the device and the warranty will be voided.

WARNING

Risk of injury: The use of parts not approved by the manufacturer can lead to personal injury and damage or malfunction of the meter or equipment. The replacement parts in this section are approved by the manufacturer.

Description	REF
Printer paper, pack of five rolls	93065
Control strip*	913470
Barcode scanner	93075
Case for QUANTOFIX® Relax	930889
Manual	930933
Mains adapter	930995
Strip holder	930890
* see Chapter 14.3 "How to run a control strip measurement"	

- 6 AA type batteries (optional)
- RS232 to LAN converter for connection to LIMS (Please contact MACHEREY-NAGEL for further information and instructions)

19. Manufacturer information

MACHEREY-NAGEL GmbH & Co. KG

If you have questions or need technical assistance after reading the manual, contact:

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The screenshot shows a screen titled "ID:" with the number "123456789" displayed. Below the display is a numeric keypad with buttons for digits 1-9, 0, and a back arrow. At the bottom, there are two large buttons: a red "X" on the left and a green checkmark on the right.

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The screenshot shows a screen titled "ID:" with the text "ABCD" displayed. Below the display is a character keypad with buttons for letters A-Z, numbers 0-9, and a back arrow. At the bottom, there are two large buttons: a red "X" on the left and a green checkmark on the right.

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