

Test strips. Reimagined.

QuantifyGo[®]

The smartphone upgrade
for QUANTOFIX[®]

Dip.
Scan.
Done!



MACHEREY-NAGEL

www.mn-net.com



Easy handling

How it works

With *QuantifyGo*[®], you can now evaluate your QUANTOFIX[®] test strips digitally and easily – directly with your smartphone and without installing an app.



- 1 Scan the QR code on the folding box or on the inside of this flyer.



- 2 Press the "Scan" button in the *QuantifyGo*[®] application.



- 3 Place the test strip (after dipping it into the solution) onto the scanning template and take a photo.

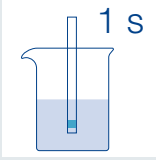
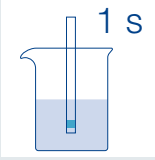
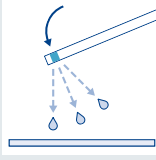
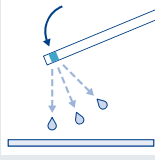


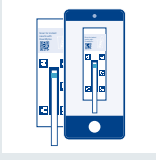
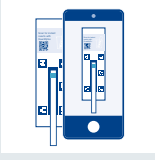


- 4 Your result appears immediately on the screen.

Test it yourself!

Experience the live digital evaluation of your QUANTOFIX[®] test strips with *QuantifyGo*[®].

Place the test strip on the designated side of the folding box or directly onto the scanning area on the inside of the flyer!

QUANTOFIX [®] Peroxide 25	QUANTOFIX [®] Peroxide100
 <p>1 s</p>	 <p>1 s</p>
	
 <p>15 s</p>	 <p>5 s</p>
	

* Note: To detect hydroperoxides in organic solvents, moisten the test area with a drop of water after the solvent has evaporated or dried.

Try it now!

Clear readings – objective, seamless, and free of charge!

- Scan the QR code with your smartphone
- Place the strip on the box graphic below
- Get your digital result instantly

Good to know

Optional registration for team features, data storage, and downloads.

Scan for instant results with QuantifyGo



Scan the QR code to go directly to QuantifyGo®!



MACHEREY-NAGEL MN
www.mn-net.com

QuantifyGo®

One tool. All data. Endless possibilities.

Instant reporting



Objective values in mg/L. Stop comparing color scales by eye.

Your data, your structure

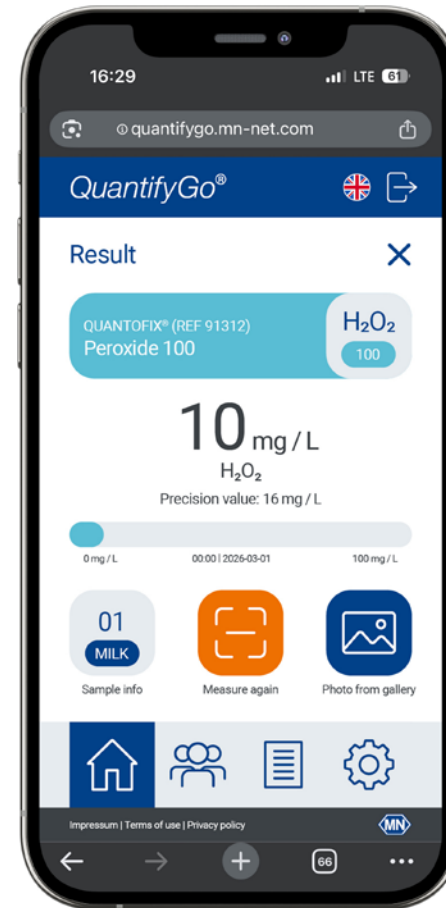


Link results directly to sample types (e.g., milk) or specific locations.

Collaborative teamwork



Share results with your team in real-time—ideal for shift handovers.



Documentation made easy



Filter measurements and export them directly as PDF or Excel reports for your audits.

Scan and go



No download, no IT approval required. Simply scan the QR code and analyze directly in your browser.

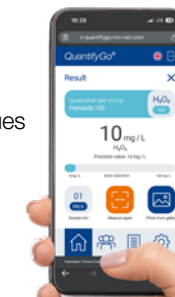
Free digital upgrade



For QUANTOFIX® users, all features—from scanning to export—are completely free of charge.

QuantifyGo® – Flexible and mobile

- Mobile analysis on-site for flexible results right where you need them.
- Digital objectivity guarantees precise values by eliminating manual reading errors.
- Digital data management enables easy archiving and seamless data export.
- An instant start without hardware allows for immediate use with no extra costs.



Ready for your digital analytics

Register to discover even more

- Save and download your results
- Add sample information
- Collaborate in teams and share data

Free!

Digitally evaluable test strips



QUANTOFIX® Peroxide 25
Measuring range 0–25 mg/L
REF 91319

QUANTOFIX® Peroxide 100
Measuring range 0–100 mg/L
REF 91312

More test strips coming soon!

Everything at a glance

Visit us online for more information, tutorials, and to download your free scanning template. Learn all about the benefits of QuantifyGo® and how to optimize your workflow.



Help shape the future!

We are developing QuantifyGo® specifically for you. Which test strip would you like to evaluate digitally next?

Submit your requested parameter:
support@mn-net.com



Your opinion matters!

Help us make the application even better. Every piece of feedback goes directly into our development process.

Any questions?

Our service team is happy to assist you:

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
Customer Service Center

Tel.: +49 24 21 969-333

support@mn-net.com



Your contact person:



MACHEREY-NAGEL GmbH & Co. KG

Valencienner Str. 11

52355 Düren · Germany