

Fluconazole and Related Substances – USP

MACHERY NAGEL application department · Dr. H. R. Wollseifen, T. Kretschmer, R. Nuessgen

Application benefits

- HPLC method with faster separations within allowable adjustments
- Shorter runtimes
- Lower solvent consumption
- Optimized system suitability

MN products

REF 760083.46

EC HPLC column (analytical),
NUCLEODUR® C18 Gravity, 3 µm,
150 × 4.6 mm

REF 763136.46

EC HPLC column (analytical),
NUCLEOSHELL® RP 18, 2.7 µm,
150 × 4.6 mm

REF 763134.46

EC HPLC column (analytical),
NUCLEOSHELL® RP 18, 2.7 µm,
100 × 4.6 mm

REF 702107

Screw closure, N 9, PP, yellow, center
hole, Silicone white/PTFE red, 1.0 mm

REF 702079

Screw neck vial, N 9, 11.6x32.0 mm,
1.5 mL, label, flat bottom, amber,
silanized

MN application numbers

HPLC: 129330

HPLC: 129340

HPLC: 129350

Keywords

Fluconazole, USP, NUCLEOSHELL®
RP18, NUCLEODUR® C18 Gravity,
L1, United States Pharmacopeia

Introduction

The USP monograph describes the separation of Fluconazole from impurities. This work starts using a fully porous HPLC phase and shows the benefits using superficially porous particles. The method optimization was performed to achieve shorter run time and system suitability results within allowable adjustments.

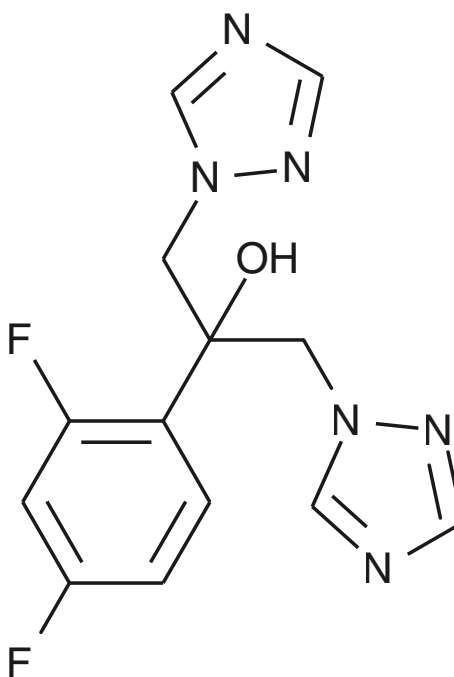


Figure 1: Structure of fluconazole.

USP Monograph: Fluconazole Method Details

Method Parameter	Description
Standard solution	10 µg/mL each of USP Fluconazole RS, USP Fluconazole Related Compound A RS, USP Fluconazole Related Compound B RS, and USP Fluconazole Related Compound C RS, dissolved in acetonitrile, and then diluted in Mobile phase.
Column size	150 × 4.6 mm
Stationary Phase	3.5-µm packing L1
Mobile Phase	Acetonitrile and water (20:80)
Flow Rate	0.5 mL/min
Temperature	40 °C
Detection	260 nm
Injection	20 µL
Run Time	1.5 times the retention time of Fluconazole (about 6 min)
Elution Order	1. Fluconazole related compound A 2. Fluconazole related compound B 3. Fluconazole related compound C 4. Fluconazole
Suitability requirements	
Resolution:	NLT 1.5 between fluconazole related compound B and fluconazole related compound C.
Relative standard deviation:	NMT 5.0 % for each peak.

* Fluconazole (USP-1271700), Fluconazole Related Compound A (USP-1271711), Fluconazole Related Compound B (USP-1271722), and Fluconazole Related Compound C (USP-1271733), were purchased from Labmix24 GmbH; Postal address: Industriestrasse 18A - 46499 Hamminkeln (Germany).

Table 1a: USP Monograph: Fluconazole Method Details

Chromatographic methodology improvements

Figure 2: a

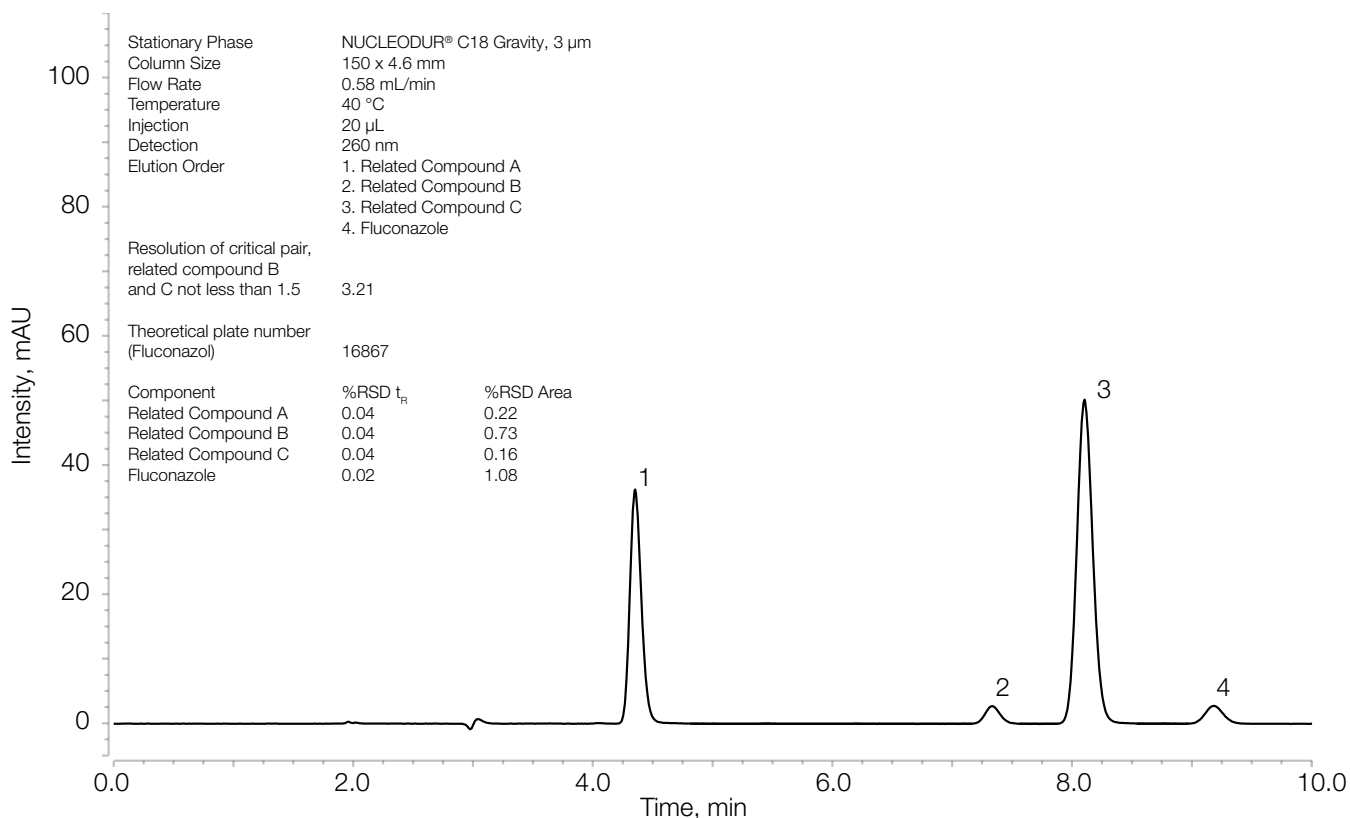


Figure 2: b

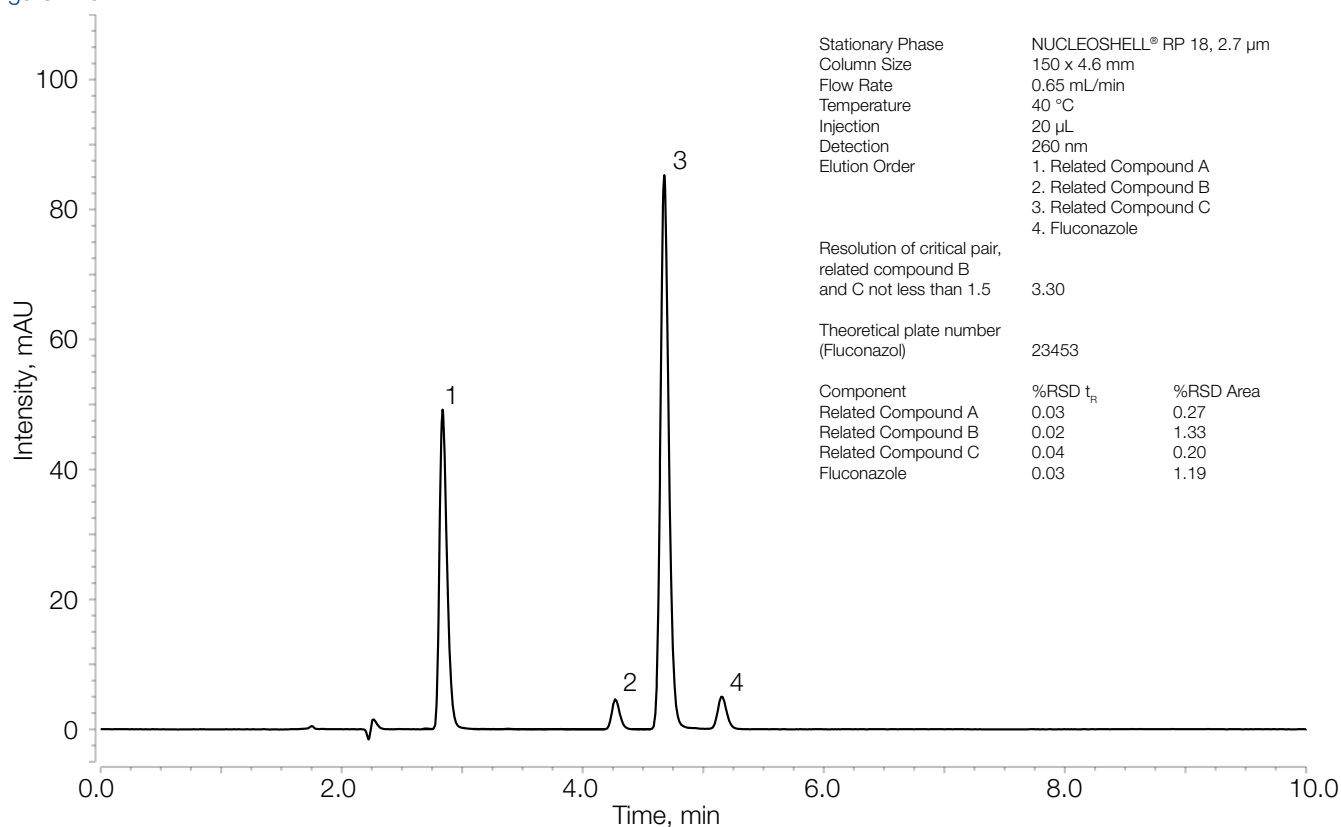


Figure 2: c

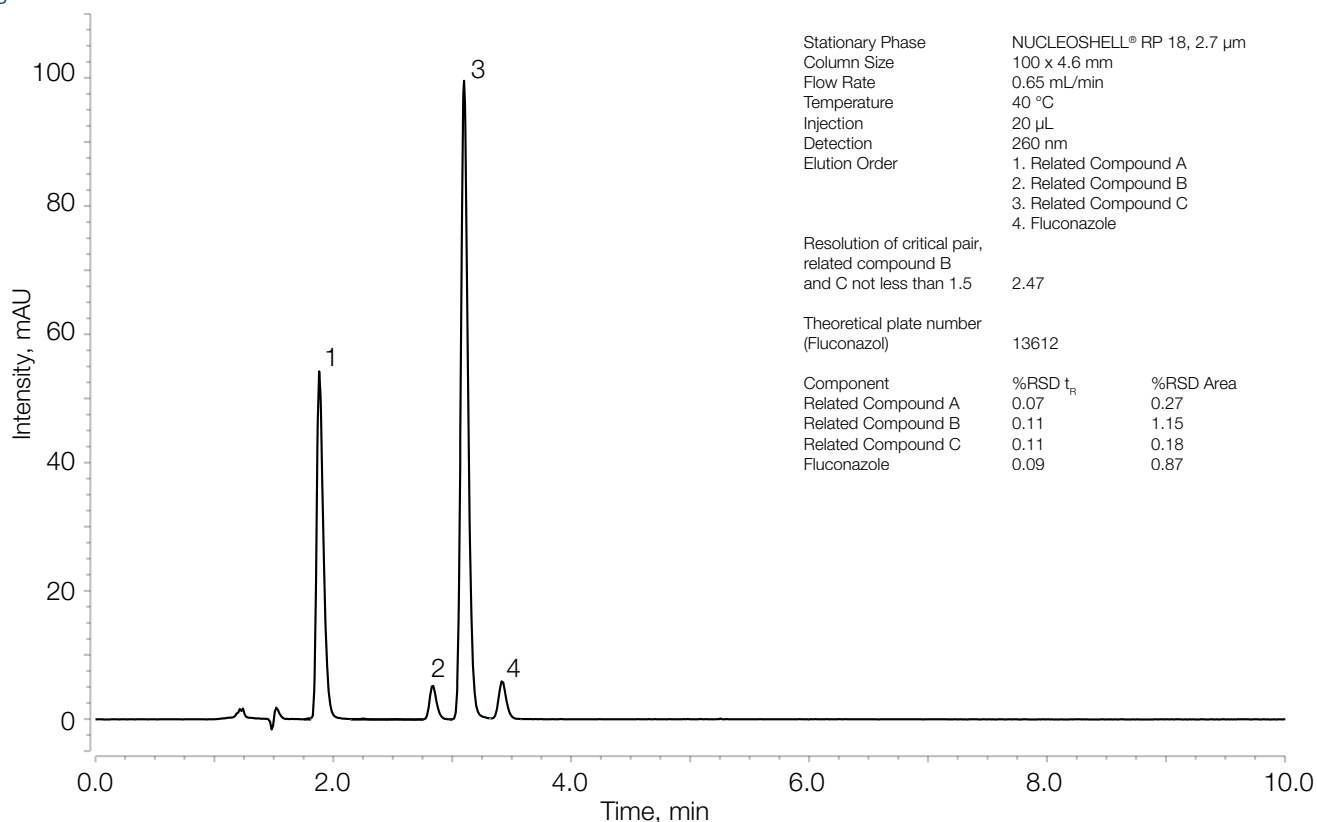


Figure 2: a: EC HPLC column (analytical), NUCLEODUR® C18 Gravity, 3 µm, 150 × 4.6 mm, b: EC HPLC column (analytical), NUCLEOSHELL® RP 18, 2.7 µm, 150 × 4.6 mm, c: EC HPLC column (analytical), NUCLEOSHELL® RP 18, 2.7 µm, 100 × 4.6 mm.

Results

Method Parameter	Allowed Adjustments (isocratic elution)*	Method 1 (figure 2a)	Method 2 (figure 2b)	Method 3 (figure 2c)
Mobile phase pH	± 0.2 units	As specified	As specified	As specified
Concentration of salts in buffer	± 10%	As specified	As specified	As specified
Composition of the mobile phase	± 30% relative; cannot exceed ± 10% absolute change; cannot be reduced to zero	As specified	As specified	As specified
Stationary phase	No change of C18 allowed	NUCLEODUR® C18 Gravity	NUCLEOSHELL® RP 18	NUCLEOSHELL® RP 18
Ratio column length/ particle size	Column length to particle size diameter ratio can be adjusted between – 25% and + 50%	150 mm / 3 µm (+ 16.7%*)	150 mm / 2.7 µm (+ 29.6%*)	100 mm / 2.7 µm (– 13.6%*)
Column internal diameter	Can be adjusted so long as linear velocity is maintained	4.6 mm as specified	4.6 mm as specified	4.6 mm as specified
Flow rate	± 50% after adjustment due to a change in column dimensions	0.58 mL/min (± 0% after adjustments)	0.65 mL/min (± 0% after adjustments)	0.65 mL/min (± 0% after adjustments)
Column temperature	± 10 °C	40 °C as specified	40 °C as specified	40 °C as specified
Injection volume	Can be adjusted as much as needed; must be consistent with linearity, precision, and detection requirements	20 µL as specified	20 µL as specified	20 µL as specified
Detection [nm]	No change permitted	260 nm as specified	260 nm as specified	260 nm as specified
Retention time Fluconazole [min]		8.101 min	5.152 min (– 36.4%**)	3.418 min (– 57.8%**)
Theoretical plate number (Fluconazol)	Within -25% to 50%, relative to the prescribed column***	16867	23453 (+ 39.1%**)	13612 (– 19.3%**)
Suitability requirements				
Resolution:	NLT 1.5 between flucon-azole related compound B and fluconazole related compound C	3.21	3.30	2.47
%RSD t _r :	NMT 5.0% for each peak.	0.02 – 0.04 (see Figure 2a)	0.02 – 0.04 (see Figure 2b)	0.07 – 0.11 (see Figure 2c)
%RSD Area:	NMT 5.0% for each peak.	0.16 – 1.08 (see Figure 2a)	0.20 – 1.33 (see Figure 2b)	0.18 – 1.15 (see Figure 2c)

* change in comparison to USP method ** change in comparison to method 1 *** column used in method 1.

Conclusion

The fully porous NUCLEODUR® C18 Gravity, 3 µm, 150 × 4.6 mm HPLC column from MACHERY NAGEL fulfills all requirements of the USP monograph (Fluconazole and Related Substances). By using superficially porous NUCLEOSHELL® analytical columns the runtime of the method can be reduced by up to 57.8% (with NUCLEOSHELL® RP 18, 2.7 µm, 100 × 4.6 mm) compared to fully porous NUCLEODUR® silica gel, while keeping all method parameters well within the allowed adjustment range of the USP monograph. The reduction in

runtime leads to a lower solvent consumption optimizing the analysis of Fluconazol with regard to the guidelines of green chemistry. We were also able to improve the peak intensity with NUCLEOSHELL® columns compared to the original method with fully porous silica gel.

MACHERY NAGEL shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material. Information, descriptions, and specifications in this publication are subject to change without notice.

www.mn-net.com

MACHERY-NAGEL



MACHERY-NAGEL GmbH & Co. KG · Valencienner Str. 11 · 52355 Düren · Germany

DE +49 24 21 969-0 info@mn-net.com

CH +41 62 388 55 00 sales-ch@mn-net.com

FR +33 388 68 22 68 sales-fr@mn-net.com

US +1 888 321 62 24 sales-us@mn-net.com