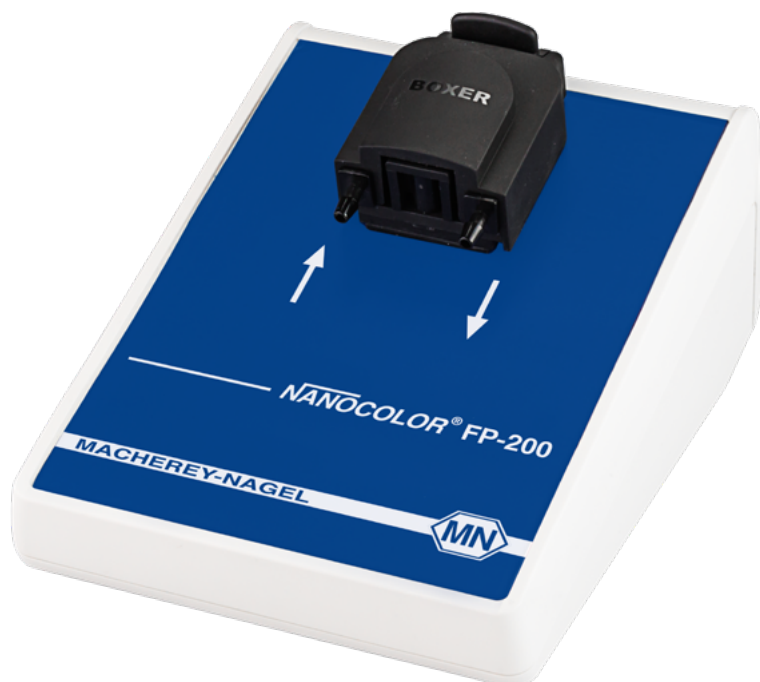
MACHEREY-NAGEL

NANOCOLOR® FP-200

- Handbuch
- Manual
- Manuel

<b>Deutsch</b> .....	<b>3</b>
<b>English</b> .....	<b>22</b>
<b>Français</b> .....	<b>41</b>
<b>Español</b> .....	<b>60</b>

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<b>1. Introduction</b> .....	<b>23</b>
<b>2. CE mark</b> .....	<b>23</b>
<b>3. Nameplate</b> .....	<b>23</b>
<b>4. Explanation of symbols</b> .....	<b>24</b>
<b>5. Safety instructions</b> .....	<b>24</b>
<b>6. Chemical and biological safety</b> .....	<b>25</b>
<b>7. Technical data</b> .....	<b>25</b>
<b>8. Setup and connection of the device</b> .....	<b>26</b>
8.1 Scope of delivery .....	26
8.2 Setting up the device.....	26
<b>9. Installation with NANOCOLOR® VIS II and <sup>UV</sup>I<sub>VIS</sub> II</b> .....	<b>27</b>
<b>10. Operation and menus</b> .....	<b>28</b>
10.1 Activating the sipper module on the spectrophotometer.....	28
10.2 Setting the pumping, rinsing, and settling times .....	28
10.3 Programming of method specific sipper module settings.....	30
10.4 Activating and deactivating the sipper module.....	31
10.5 Executing a stored program .....	32
10.6 Taking a single wavelength or multi wavelength measurement.....	33
10.7 Performing a scan .....	34
10.8 Performing a color measurement.....	35
<b>11. Maintenance and cleanup of the device</b> .....	<b>36</b>
11.1 Cleaning the flow cuvettes .....	36
11.2 Cleaning the hoses .....	36
11.3 Replacing the pump hose .....	37
<b>12. Errors, causes, and solutions</b> .....	<b>38</b>
<b>13. Service</b> .....	<b>39</b>
13.1 Replacement parts and accessories.....	39
13.2 Disposal .....	39
13.3 Warranty.....	40
13.4 Contact.....	40

# 1. Introduction

The *NANOCOLOR*<sup>®</sup> FP-200 is an optional accessory for the *NANOCOLOR*<sup>®</sup> VIS II and *NANOCOLOR*<sup>®</sup> UV/VIS II spectrophotometers. Using a peristaltic pump, a constant amount of liquid is pumped through a flow cuvette and the measurement is started automatically on the photometer. It is possible to improve the measurement accuracy by working with the *NANOCOLOR*<sup>®</sup> FP-200. The same optical conditions apply to the zero measurement and the measurement of the sample. Errors due to optical differences between different cuvettes are avoided; all solutions to be analyzed flow through one cuvette.

# 2. CE mark

The CE mark declares that the product complies with the harmonization legislation of the European Community listed below:

- European Directive 2011/65/EU on the restriction of the use of certain hazardous substances in electrical and electronic equipment (RoHS 2)
- European Directive 2012/19/EU on waste electrical and electronic equipment (WEEE)
- European Directive 2014/30/EU on the harmonization of the laws of the member states relating to electromagnetic compatibility (EMC)

# 3. Nameplate

## 4. Explanation of symbols

The nameplate of the device contains the symbols and terms listed below with the following meanings:

Term / Symbol	Meaning
NANOCOLOR® FP-200	Device type designation
SN	Serial number of the device
5 V, 500 mA	Power supply
	Warning against biological hazards to users that can result from the use of the device. Note also the information in section 6 “Chemical and biological safety”.
	Caution (note the information in the documentation)! Follow the safety instructions in the operating manual of the device.
	According to 2012/19/EU, it is prohibited to dispose of the device through public waste disposal systems. Note also the information in section 13.2 “Disposal”.
	The CE symbol indicates fulfillment of the applicable harmonization legislation of the European Community.
MACHEREY-NAGEL 	Identification of the manufacturer

## 5. Safety instructions

*Caution:* Follow the instructions in this section carefully to ensure the device does not become damaged.

**Warning:** The instructions in this section must be followed carefully to ensure that no further dangers or hazards exist, otherwise the device will very likely become damaged.

Note: Information, which illustrate aspects from the main text.

*Caution:* The NANOCOLOR® FP-200 may only be used as specified by the manufacturer and only be correspondingly qualified personnel. Note all markings and signs mounted on the device or stated in the operating manual. Ignoring them can lead to personal injuries and/or damage to the device.

*Caution:* The NANOCOLOR® FP-200 is delivered with a plastic hose adapters. The plastic hose adapter is only suitable for use of MACHEREY-NAGEL rectangular cuvette tests. Resistance to aggressive chemicals is not guaranteed. Please use a stainless steel hose adapter in this case.

## 6. Chemical and biological safety

**Warning:** Working with chemicals can be hazardous. Handling chemicals without protection can lead to serious injuries. When working with the *NANOCOLOR*<sup>®</sup> FP-200, and especially when removing the hoses, wear the required personal protective equipment.

During normal operation of this device, it may be necessary to use chemicals that are hazardous to your health or biologically harmful samples.

Before handling these substances, read all hazard statements and safety information printed on the containers with the original solutions and in the safety data sheets. All used solutions must be disposed of in accordance with the national laws and regulations. The type of protective equipment must be selected according to the concentration and amount of the hazardous substance present at the corresponding workplace. Rinse the hoses and cuvettes thoroughly after use with distilled water or a rinsing solution recommended by the manufacturer.

## 7. Technical data

Specification	Description
Type	Peristaltic pump
Operation	Fully automatically using the <i>NANOCOLOR</i> <sup>®</sup> spectrophotometer
Wavelength range	Plastic: 340 nm–900 nm Quartz glass: 190 nm–900 nm
Power supply	Via USB cable, 5 V, 500 mA
Rinsing volume	min. 10 mL
Flow rate	1 mL/s
Interface	USB B
Ambient conditions (transport, storage)	-10 °C–60 °C, 20%–80% relative humidity (without condensate formation)
Ambient conditions (operation)	10 °C–40 °C, 20%–80% relative humidity (without condensate formation)
Dimensions	125 mm x 177 mm x 69 mm
Weight	365 g
Enclosure protection class	IP30
Certification	CE-certified

## 8. Setup and connection of the device

### 8.1 Scope of delivery

Check the delivery for completeness. If some of the scope of delivery is missing, then contact MACHEREY-NAGEL directly or your sales partner.

- Sipper module *NANOCOLOR*<sup>®</sup> FP-200, pre-mounted pump hose
- USB connecting cable
- Tygon<sup>®</sup> hose, 2 m, ID = 3 mm
- Tygon<sup>®</sup> hose, 2 m, ID = 1.6 mm
- 8 reducers
- Plastic hose adapter
- Operating manual (these instructions)

### 8.2 Setting up the device

*Caution:* Place the device on a solid, level surface. Do not move the device while taking measurements. Set up the device in a work area where there is little dust.

*Caution:* When setting up the *NANOCOLOR*<sup>®</sup> FP-200, make sure that the hoses are not kinked. Always check to ensure the waste container is not full.

*Caution:* Please use the USB connecting cable provided to connect the device. You cannot use a USB hub for connection.

*Caution:* There can always be residual liquid in the hoses. When working with the *NANOCOLOR*<sup>®</sup> FP-200, and especially when removing the hoses, wear the required personal protective equipment.

## 9. Installation with **NANOCOLOR® VIS II** and **UV/VIS II**

1. The pump head is not fastened in place yet upon delivery of the **NANOCOLOR® FP-200**. Flip the lever on the back end of the peristaltic pump towards the front before using the device.
2. Place the **NANOCOLOR® FP-200** to the right of the photometer.
3. Connect the sipper module to one of the two USB A interfaces on the photometer using the connecting cable provided.
4. Remove the cover from the cuvette slot of the **NANOCOLOR® UV/VIS II** or **NANOCOLOR® VIS II**.
5. Place the desired flow cuvette in the cuvette slot of the photometer. Clean the cuvette thoroughly with a lint-free cloth before inserting it in the slot.
6. Cut a max. 40 cm long piece from the hose provided and use it to connect the hose adapter to the inlet nozzle of the flow cuvette.
7. Cut a max. 40 cm long piece from the hose provided and use it to connect the outlet nozzle of the flow cuvette to the input of the peristaltic pump.
8. Connect a part of the remaining hose to the output of the peristaltic pump and place the other end in a suitable waste container.
9. Activate the sipper module on the spectrophotometer (section 10.1)

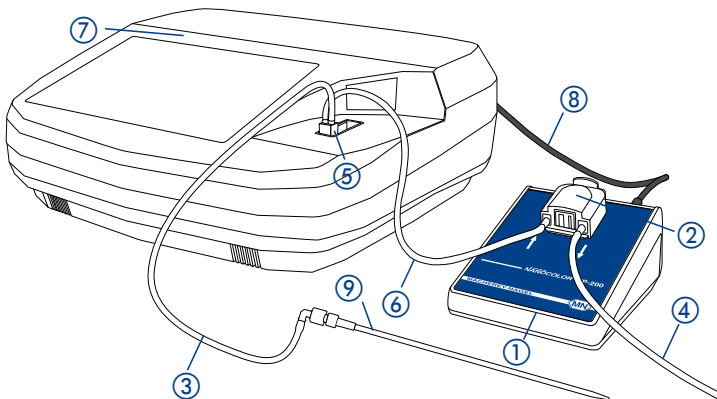
*Caution:* When taking measurements in the wavelength range of 190 nm–340 nm, place the special cuvette slot cover on the **NANOCOLOR® UV/VIS II** (REF 919610).

*Caution:* Make sure that the hoses are not pinched. The pump volume of the setup can be optimized by shortening the hoses.

*Caution:* The thin hoses are not suited for the direct connection to cuvettes or the Sipper module. Please use the reducers and a short part of the thick hose for connection.

- ① Sipper module *NANOCOLOR*<sup>®</sup> FP-200
- ② Peristaltic pump
- ③ Supply hose
- ④ Drain hose
- ⑤ Flow cuvette
- ⑥ Connecting hose
- ⑦ Photometer
- ⑧ USB connecting cable
- ⑨ Hose adapter

## 10. Operation and menus

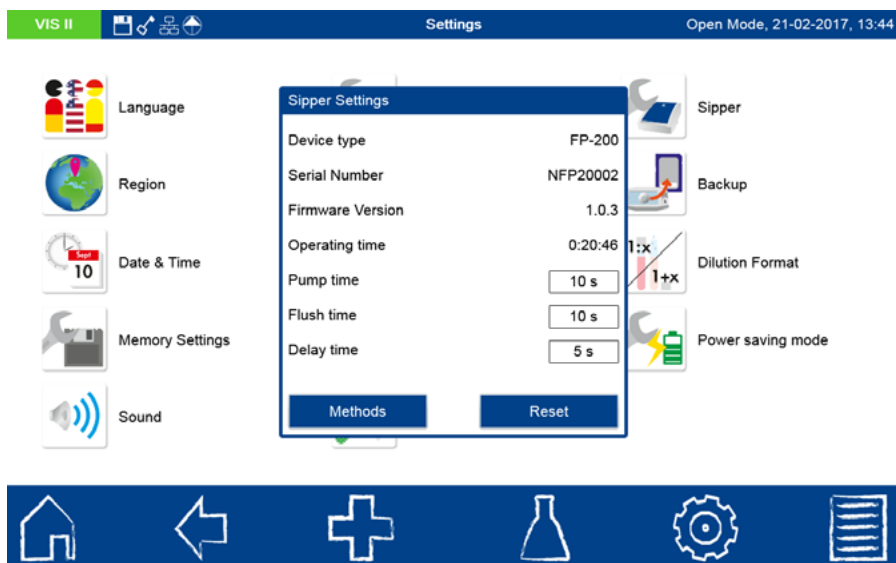
### 10.1 Activating the sipper module on the spectrophotometer

After connecting the sipper module to the spectrophotometer, open the connectivity menu via  →  → . Activate the checkbox for the FP-200 and the USB A interface. Please confirm the changes by pressing . The sipper module is detected automatically by the spectrophotometer. The  symbol appears in the status bar of the photometer.

### 10.2 Setting the pumping, rinsing, and settling times

The sipper module settings are configured in a separate settings menu on the spectrophotometer. After connecting the sipper module to the spectrophotometer, open the configuration menu via  →  → .

A window displaying the properties of the sipper module appears.

Pumping time:

The pumping time determines the sample volume that will be pumped through the flow cuvette. To ensure a 50 mm flow cuvette is completely rinsed, the pumping time must be set to a min. of 10 s.

*Caution:* The pumping time should be determined individually for each flow cuvette and each measurement setup. To do this, you can measure how an air bubble wanders through the hose, for example, or measure the absorbance continuously for two solutions with different concentrations. Pumping time, rinsing time and settling time can be set method specific (section 10.3).

- Rinsing time: The rinsing time determines the volume of rinsing solution that will be pumped through the flow cuvette after every measurement. Rinsing can be performed manually, if necessary after each measurement.
- Settling time: The settling time determines the wait time between the pumping operation and the measurement. The settling time allows air bubbles that have formed during the pumping operation to escape and allows time for turbulence in the sample to stop.
- Serial number: Shows the serial number of the connected sipper module.
- Device type: Shows the type of device connected.
- Firmware version: Shows the firmware version of the connected sipper module.
- Hours of operation: Shows the number of hours of operation of the connected sipper module.

The times are entered in seconds using the numeric keypad. The keypad opens by clicking on the corresponding entry.

The „Reset“ function enables a reset of the pumping time, rinsing time and settling time to factory settings.


The field “Methods” will open the menu for programming the sipper methods (section 10.3).

### 10.3 Programming of method specific sipper module settings

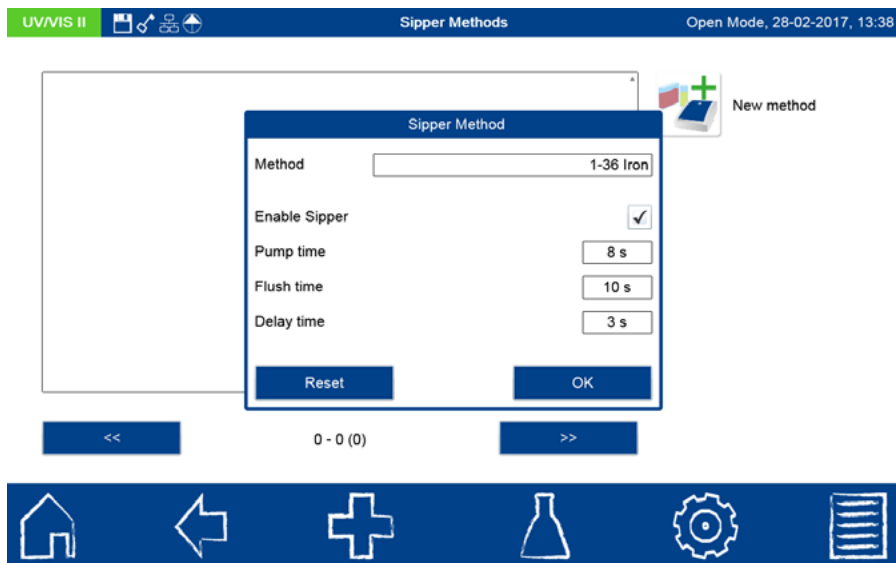Before using the sipper module the methods to be run with the sipper module must be defined. This allows for the use of different parameters for pumping time, rinsing time and settling time for each method. These settings are stored method specific by the spectrophotometer.

After connection of the sipper module to the spectrophotometer, please open the configuration menu via  →  → .

The popup with the sipper settings will appear (section 10.2). Chose the field “Methods” to enter the menu for programming the method specific sipper module settings. Clicking on the icon  will open a popup window for creation of a new sipper method.

Click on the field “Method” to choose the desired method for the programming. A list with all methods suitable for evaluation in rectangular cuvettes will open. Please choose the method of choice from the list. Set the pumping time, the rinsing time and the settling time. The times are entered in seconds using the numeric keypad. The keypad opens by clicking on the corresponding entry.

The function „Reset“ will reset the settings for pumping time, rinsing time and settling time to the values set in the general sipper settings (section 10.2).

Activating the checkbox “Enable Sipper” will automatically activate the sipper module when calling up the respective method. Confirming with  stores the new method in the list of the sipper methods.

After marking an entry from the list of sipper methods the options on the right hand side become available. Confirming  will open the menu for setting the method specific sipper settings. The settings of the method can now be edited.

Confirming the icon  will delete the selected method after confirming a safety question.

The checkbox behind the method will activate or deactivate the automatic use of the sipper when calling up the respective method.

When calling up a method for performing a measurement, the spectrophotometer will detect automatically, if the method is programmed for the use with the sipper module. The method specific sipper module settings will be called up automatically.

Note: If no methods are defined in the list of the sipper methods, the sipper module will not be used even when connected and activated.

#### 10.4 Activating and deactivating the sipper module

In the menus for scan and color measurement, the sipper module can be deactivated via the measurement menu to enable the measurement of samples without the sipper module even when connected and activated. To deactivate the sipper module, press  in the corresponding measurement menu and select “Deactivate sipper”. To reactivate the sipper module, proceed in the same manner and then select “Activate sipper”.

## 10.5 Executing a stored program

Call the desired test using its test number or from the test selection list. When the sipper module is activated, you will see the  icon in the status bar of the photometer. In case methods are programmed in the list of sipper methods (section 10.3), the settings stored with the method will be automatically called used.

Note: Please consider the information in the instructions for use of the respective test kit.

Hold the intake hose in the zero solution and press the  icon. The zero solution is pumped into the flow cuvette and the zero measurement is taken automatically after the settling time has expired. The time remaining until the measurement will be taken is shown in the display. You can cancel the measurement by pressing .

Hold the intake hose in the sample solution to be analyzed and press the  icon. The sample solution is pumped into the flow cuvette and the sample measurement is taken automatically after the settling time has expired. The time remaining until the measurement will be taken is shown in the display. You can cancel the measurement by pressing .

The flow cuvette is rinsed using distilled water or the next sample solution. The rinsing operation is initiated in the corresponding measurement menu by pressing . The result of the measurement is displayed in the result window.

1161 Chlor

4,0 mg/L Cl<sub>2</sub>

540 nm IQC

4,0 mg/L Cl<sub>2</sub>

0,737 A  
18,3% T

Sample Type Sample Author Location Comment Flush Zero Measure

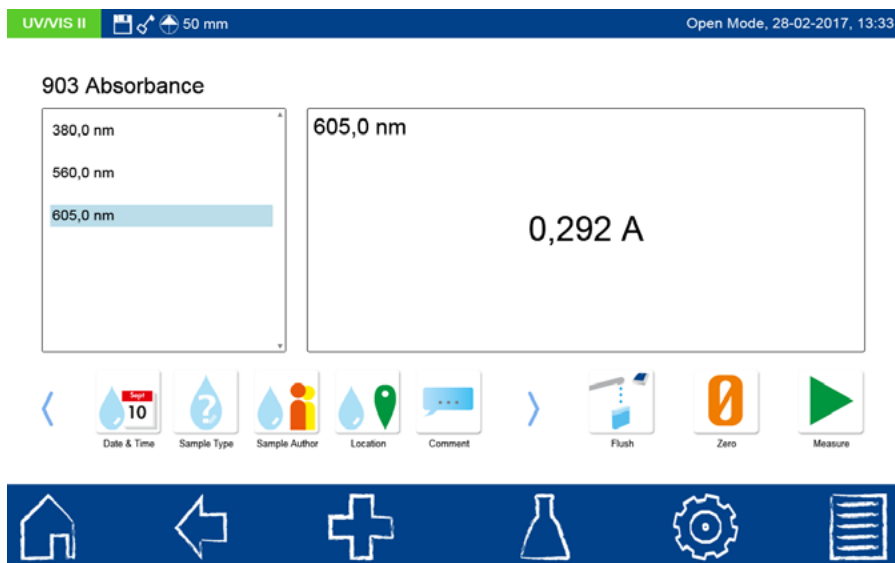
You can enter various information on the sample using the icons below the result window. The information entered is saved when you exit the menu or when the next measurement is taken.

## 10.6 Taking a single wavelength or multi wavelength measurement

Open the desired menu via  → . Enter the required information and confirm with . Hold the intake hose in the zero solution and press the  icon. The zero solution is pumped into the flow cuvette and the zero measurement is taken automatically after the settling time has expired. The time remaining until the measurement will be taken is shown in the display. You can cancel the measurement by pressing .

Hold the intake hose in the sample solution to be analyzed and press the  icon. The sample solution is pumped into the flow cuvette and the sample measurement is taken automatically after the settling time has expired. The time remaining until the measurement will be taken is shown in the display. You can cancel the measurement by pressing .

The flow cuvette is rinsed using distilled water or the next sample solution. The rinsing operation is initiated in the corresponding measurement menu by pressing . The result of the measurement is displayed in the result window.

The screenshot displays the instrument's interface. At the top, a status bar shows "UV/VIS II" on the left and "Open Mode, 28-02-2017, 13:33" on the right. Below this, the main display area is titled "903 Absorbance". On the left side of the display, there is a list of wavelengths: 380,0 nm, 560,0 nm, and 605,0 nm. The 605,0 nm option is highlighted with a blue bar. To the right of this list, the selected wavelength "605,0 nm" is displayed. In the center of the display, the absorbance value "0,292 A" is shown. Below the main display area, there is a row of nine icons for data entry: a left arrow, "Date & Time" (with a calendar icon and "10"), "Sample Type" (with a question mark icon), "Sample Author" (with a person icon), "Location" (with a location pin icon), "Comment" (with a speech bubble icon), a right arrow, "Flush" (with a flask icon), "Zero" (with a zero icon), and "Measure" (with a play icon). At the bottom of the screen, there is a dark blue navigation bar with six white icons: a home icon, a back arrow, a plus sign, a flask icon, a gear icon, and a list icon.

You can enter various information on the sample using the icons below the result window. The information entered is saved when you exit the menu or when the next measurement is taken.

## 10.7 Performing a scan

Open the menu for scans via  → . Enter the required information and confirm with . Hold the intake hose in the zero solution and press the  icon. The zero solution is pumped into the flow cuvette and the zero measurement is taken automatically after the settling time has expired. The time remaining until the measurement will be taken is shown in the display. You can cancel the measurement by pressing .

Hold the intake hose in the sample solution to be analyzed and press the  icon. The sample solution is pumped into the flow cuvette and the sample measurement is taken automatically after the settling time has expired. The time remaining until the measurement will be taken is shown in the display. You can cancel the measurement by pressing .

The flow cuvette is rinsed using distilled water or the next sample solution. The rinsing operation is initiated in the corresponding measurement menu by pressing . After the scan is complete, the results are displayed as a curve in the result window.

You can enter various information on the sample using the icons below the result window. The information entered is saved when you exit the menu or when the next measurement is taken. The scan can be edited using the scan analysis and the  icon (see the NANOCOLOR® VIS II and <sup>UV</sup>/VIS II spectrophotometer manual).

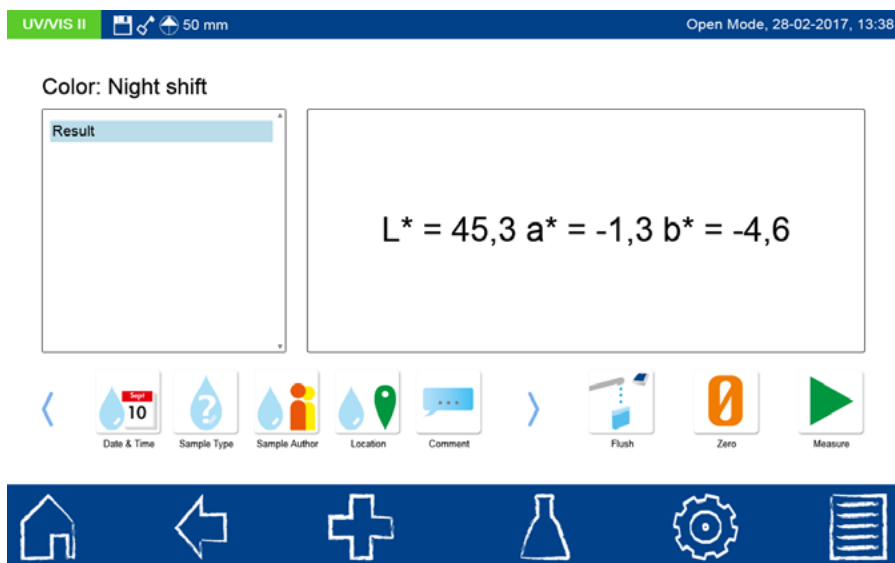
## 10.8 Performing a color measurement

Open the menu for color measurements using  → . Select the desired color measurement. Enter the required information and confirm with .

Hold the intake hose in the zero solution and press the  icon. The zero solution is pumped into the flow cuvette and the zero measurement is taken automatically after the settling time has expired. The time remaining until the measurement will be taken is shown in the display. You can cancel the measurement by pressing .

Hold the intake hose in the sample solution to be analyzed and press the  icon. The sample solution is pumped into the flow cuvette and the sample measurement is taken automatically after the settling time has expired. The time remaining until the measurement will be taken is shown in the display. You can cancel the measurement by pressing .

The flow cuvette is rinsed using distilled water or the next sample solution. The rinsing operation is initiated in the corresponding measurement menu by pressing . After the color measurement has been taken, the result is shown in the result window.

The screenshot shows the instrument's display in 'Open Mode' on 28-02-2017 at 13:38. The display is titled 'Color: Night shift'. On the left, there is a 'Result' window. On the right, the color measurement results are displayed:  $L^* = 45,3$ ,  $a^* = -1,3$ , and  $b^* = -4,6$ . Below the result window is a row of icons for sample information: Date & Time (showing '10'), Sample Type (question mark), Sample Author (person icon), Location (map pin), Comment (speech bubble), Flush (water tap), Zero (orange circle with '0'), and Measure (green play button). At the bottom of the screen is a dark blue navigation bar with icons for Home, Back, Add, Flask, Settings, and List.

You can enter various information on the sample information using the icons below the Result window. The information entered is saved when you exit the menu or when the next measurement is taken. The color measurement can be edited using the color analysis and the  icon (see the *NANOCOLOR<sup>®</sup> VIS II* and *UV<sub>VIS</sub> II* spectrophotometer manual).

## 11. Maintenance and cleanup of the device

*Caution:* The work described in this chapter may only be performed by qualified personnel.

*Caution:* The housing of the device can be cleaned with a moist cloth. Immediately wipe off all liquids splashed on the device. Do not use solvents to clean the device.

### 11.1 Cleaning the flow cuvettes

The flow cuvettes should be rinsed with distilled water at the beginning and the end of every series of measurements. If they are very dirty, then they should be rinsed out several times. Empty the flow cuvettes, if not used or stored for a longer period of time by feeding air through the cuvette. Hold the cuvette with the outlet nozzle pointing downwards at an angle to empty it. If the cuvette is very dirty, then please follow the cleaning instructions from the manufacturer of the flow cuvettes.

### 11.2 Cleaning the hoses

*Caution:* There can always be residual liquid in the hoses. When working with the NANOCOLOR® FP-200, and especially when removing the hoses, wear the required personal protective equipment.

The hoses should be rinsed with distilled water at the beginning and the end of every series of measurements. If they are very dirty, then they should be rinsed out several times. Empty the hoses if not used or stored for a longer period of time by feeding air through the hoses.

*Caution:* Remove the pump head after each use (section 11.3, point 4), to avoid mechanical stress to the hoses. This is the only way to guarantee a long service life.

**Warning:** The hoses are exposed to chemicals and must be replaced every 12 months at a minimum.

### 11.3 Replacing the pump hose

The pump hose is exposed to mechanical loads and chemicals and must be replaced every 12 months at a minimum.

1. Remove all residual liquids from the system by repeating the rinsing operation with air several times.
2. Disconnect the sipper module from the photometer.
3. Disconnect the hoses from the input and output of the peristaltic pump.
4. Push the lever on the back of the peristaltic pump towards the rear. The head of the peristaltic pump moves towards the rear when the lever is pushed.
5. Lift the pump head upwards and remove the pump hose and connection nozzle. *Note: Please do not dispose of the hose and the nozzles in household waste.*
6. Insert the new pump hose in the pump head and then place the pump head on the peristaltic pump.
7. Push the lever on the back of the peristaltic pump towards the front again. The cover of the peristaltic pump moves towards the front when the lever is pushed.
8. Connect the hoses back to the input and output of the peristaltic pump.
9. Reconnect the sipper module to the photometer.

## 12. Errors, causes, and solutions

Potential errors in the *NANOCOLOR*<sup>®</sup> FP-200 are displayed on the photometer.

<b>Error</b>	<b>Cause</b>	<b>Solution</b>
Sipper module not connected! Please check the USB cable connection.	Sipper is not recognized by the spectrophotometer.	Cable is not installed correctly. Wrong cable was used.
Please check the sipper and the hoses.	Pump is blocked.	Pump head not placed correctly on pump.
Please check the pump head.	Hose is defective. Wrong hose was used. Pumping head not installed correctly.	Replace the pinched hose. Use only original hoses and adapters. Install pump head as described in the user manual.
The Sipper module does not pump liquid.	The system has a leakage and sucks air.	Please control all connections for correct fixing.
Hardware error	Sipper module defective.	Please contact MACHEREY-NAGEL or your local distributor.
The sipper module is not detected by the spectrophotometer.	Connectivity settings are not set correctly. USB cable is not connected correctly or cable is broken.	Please check the connectivity settings as described in section 10.1. Please check the USB cable connection for correct contact.
The spectrophotometer shows the icon for removal of the cuvette during the starting sequence.	The flow cell is still placed in the cuvette slot of the spectrophotometer. The device can only be started with an empty cuvette slot.	Please remove the flow cell from the before starting the spectrophotometer.
Missing response from sipper	Communication between sipper module and spectrophotometer is interrupted.	Please disconnect USB connection cable and reconnect again.
Sipper module is connected and activated, but cannot be used during measurement.	There are no sipper methods defined.	Please define one or more sipper methods via the settings menu of the sipper module (section 10.3)

In case of a persisting error, please contact MACHEREY-NAGEL or your local distributor.

## 13. Service

### 13.1 Replacement parts and accessories

Item	REF
NANOCOLOR® FP-200 for NANOCOLOR® UV/VIS II and NANOCOLOR® VIS II	919180
Support stand for NANOCOLOR® FP-100 and NANOCOLOR® FP-200	919143
Plastic hose adapter for NANOCOLOR® FP-100 and NANOCOLOR® FP-200	919142
Cuvette slot cover for NANOCOLOR® UV/VIS II for measurements with the FP-200	919610
USB connecting cable for NANOCOLOR® FP-200	919687
Set of hoses for NANOCOLOR® FP-200	919181
NANOCOLOR® FP-200 manual	919182
Flow cuvette, quartz glass, 10 mm optical path, for NANOCOLOR® UV/VIS II	919626
Flow cuvette, quartz glass, 50 mm optical path, for NANOCOLOR® UV/VIS	919129
Flow cuvette, quartz glass, 50 mm optical path, for NANOCOLOR® UV/VIS II, NANOCOLOR® VIS II, and NANOCOLOR® VIS	919149
Flow cuvette, quartz glass, 2 mm optical path, for NANOCOLOR® UV/VIS II and NANOCOLOR® VIS II	919127
Flow cuvette, quartz glass, 10 mm optical path, for NANOCOLOR® UV/VIS II, NANOCOLOR® VIS II, and NANOCOLOR® VIS	919158
Fixing for 10 mm cuvette, for NANOCOLOR® UV/VIS II, NANOCOLOR® VIS II, and NANOCOLOR® UV/VIS	919136

### 13.2 Disposal

charge.

Dispose of according to Directive 2012/19/EU (WEEE). In agreement with the European Directive 2012/19/EU on waste electrical and electronic equipment (WEEE), MACHEREY-NAGEL will take back the old equipment and dispose of it free of

Note: It is prohibited to dispose of the device through public waste disposal systems. Please contact your local MACHEREY-NAGEL representative.

### **13.3 Warranty**

The warranty for this device has a duration of 24 months from the date of purchase. The original copy of the bill serves as verification and must be submitted when filling a warranty claim. The warranty is void if the device is handled and/or maintained improperly.

The warranty is limited to the repair of defective parts or - at our sole discretion - to the delivery of an error-free replacement device. The warranty period of 24 months is not affected when a warranty claim is submitted during this period. There is no right of withdrawal. All other warranty claims are excluded. This includes in particular all claims for damages resulting from consequential damage or indirect damage. In addition, the currently valid version of our general terms of sale and delivery apply as printed on all price lists.

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MACHEREY-NAGEL GmbH & Co. KG