# NANOCOLOR Ringversuch Metalle

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# **NANOCOLOR** Interlaboratory Trial Metals

## **System for Analytical Quality Control**

This **NANOCOLOR** unit consists of three interlaboratory trial solutions, as well as a PDF evaluation sheet or online form. Both are available at [www.mn-net.com/Auswertebogen925002](http://www.mn-net.com/Auswertebogen925002).

**Shelf Life:** 3 months

The Interlaboratory Trial solutions consist of multistandards of unknown composition which are checked by a neutral and accredited control laboratory. The solution is acidified with nitric acid for preservation (pH 3–5).

The Interlaboratory Trial solutions have been developed for **internal quality control** (IQC) (according to the German DWA leaflet A 704) to check the working method and equipment as well as to control the proper handling. The measured values are examined and certificated by the Interlaboratory Trial organizer.

**Recommended frequency of application:** 1x per year

### **Contents:**

1 x 30 mL Interlaboratory Trial Metals solution A

1 x 30 mL Interlaboratory Trial Metals solution B

1 x 30 mL Interlaboratory Trial Metals solution C

for analytical control of the parameters aluminium, cadmium, chromate, total chromium, copper, iron, nickel, lead and zinc

### **Hazard warning:**

Information regarding safety can be found on the box' label and in the safety data sheet. You can download the SDS from [www.mn-net.com/SDS](http://www.mn-net.com/SDS).

### **Required accessories:**

Photometer, heating block, tube tests, piston pipettes with disposable tips

### **Procedure:**

For the selection of the parameters and the **NANOCOLOR**<sup>®</sup> tube tests, the MACHEREY-NAGEL test no. and the expected concentration range are indicated on the PDF evaluation sheet.

Use the Interlaboratory Trial solution instead of a water sample.

Perform the analyses as usual following the instructions given in the manual.

We recommend to conduct the Interlaboratory Trial solution within 14 days after receipt. The analyses should be carried out within 2 days. Each specified parameter can be determined at least twice.

Document the measurement results, the reagents used and the photometer on the PDF evaluation sheet that can be filled out digitally. Save the completed PDF document and send it by e-mail to: [CSC@mn-net.com](mailto:CSC@mn-net.com). Alternatively, please use our online form at [www.mn-net.com/Auswertebogen925002](http://www.mn-net.com/Auswertebogen925002).

Afterwards, you will receive a complete evaluation of the Interlaboratory Trial accompanied with a certification of your measurement. If necessary, individual advice for improvement of the results will be provided.

Your analysis results will of course be treated confidentially and will neither be passed on nor published.

### **NANOCOLOR Interlaboratory Trial with tests and photometers of other manufacturers:**

Use the Interlaboratory Trial solution instead of a water sample in each test. Perform the analyses as usual following the manufacturer's instructions. It may be necessary to adjust the pH value before the determination.

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