

Customer information

18.05.2021, version 1, RaKr

Operating system update for spectrophotometers NANOCOLOR® VIS II and NANOCOLOR® UV/VIS II

Overview

Photometer	NANOCOLOR® VIS II and NANOCOLOR® UV/VIS II
Version number	7.0.2864.4
Availability	from 21 st of May 2021
Price	free of charge
Required accessories	USB-stick
Other requirements	Software version ≥ 1.9.0




Important note

An update of the operating system to the version mentioned above is necessary if you want to connect the new sipper pump FP-200 or if you are planning to integrate the device in a network. We recommend a backup of all photometer data before performing the operating system update. To do this, please proceed as described in the instruments user manual.



The operating system update can lead to damage of the device if the notes and instructions for the update are not followed.

Update procedure

- Extract this ZIP file to a folder of your choice on your computer.
- Copy the .arc-update file to the upper level of an USB-stick and connect the USB stick to the device.
- Press  →  →  and choose the entry "Operating system update".






The entry "Operating system update" is only available in open mode or a user with administration rights.

- Confirm the prompt with "Yes", if you want to perform the operating system update. The update process may take several minutes.



Please do not switch off your device during the update process. Please do not remove the USB stick during the update process. The device may take damage otherwise.

- Reboot the device on prompt. Wait until the device is completely rebooted and has finished to perform its self-tests. Please do not switch off the device during this process.
- Please check, if the update process was successful via  →  → . The new operating software version will now be shown.

Modifications

- Bugfixes

Contact

If you have further questions, please do not hesitate to contact us:

Technical support and customer service	Tel.:	+49 24 21 969-332
	E-Mail:	csc@mn-net.com
For orders (international)	E-Mail:	sales@mn-net.com

www.mn-net.com



MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6-8 · 52355 Düren · Germany

DE/international:

Tel.: +49 24 21 969-0

Fax: +49 24 21 969-199

E-mail: info@mn-net.com

CH:

Tel.: +41 62 388 55 00

Fax: +41 62 388 55 05

E-mail: sales-ch@mn-net.com

FR:

Tel.: +33 388 68 22 68

Fax: +33 388 51 76 88

E-mail: sales-fr@mn-net.com

US:

Tel.: +1 484 821 0984

Fax: +1 484 821 1272

E-mail: sales-us@mn-net.com