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BioFix® Lumi „Single-Shot“ Luminescent Bacteria for individual determinations in accordance with operating control and self supervision

Method:

Determination of the acute toxicity with the aid of **freeze-dried** luminescent bacteria. The measurable variable is the natural luminescence of the test microorganism *Vibrio fischeri* NRRL B-11177. Evaluated is the inhibition of the luminous intensity in the sample in comparison with an uninhibited control. The amount of inhibition of the luminous intensity in the sample is a measure of its toxicity.

Field of application:

Waste water, aqueous extracts and leachates, fresh water (surface water and ground water), sea-water, brackish water, eluates of sediments (fresh water brackish water and seawater), interstitial pore water, and single substances dissolved in water.

Measuring range:

0 – 100 % inhibition

Indication of the results:

- **% Inhibition** of the light output in the sample in comparison to an uninhibited control
- **EC_{xx}**: concentration of a sample which causes an inhibition of the light output of exactly xx % (e.g. EC₅₀ value: concentration of the sample which causes an inhibition of the light output of 50 %):
- **TU values** („Toxicity units“ in accordance with the definition of the United States Environmental Protection Agency U.S. EPA): 100 divided by EC₅₀ value.

Content of reagent sample

20 vials with freeze-dried BioFix® Lumi luminescent bacteria

20 glass cuvettes, 50 × 12 mm

1 flask with 30 mL „BioFix® Lumi Reactivation solution for „Single-Shot“ luminescent bacteria“

1 flask with 15 mL „BioFix® Lumi Control solution“

Safety precautions:

This test kit does not contain any hazardous substances for which labelling is obligatory. The luminescent microorganism *Vibrio fischeri* NRRL B-11177 has never appeared as a pathogen up to now. In accordance with the code of practice B006 1/92 ZH 1/346 of the employer's liability insurance association of the chemical industry *Vibrio fischeri* is assigned to the risk group 1 i.e. there is no risk for humans and vertebrates.

Storage:

The vials with freeze-dried BioFix® Lumi luminescent bacteria can be stored until the expiration date indicated on the packaging at temperatures of -18 to -22 °C. „BioFix® Lumi Reactivation solution for „Single-Shot“ luminescent bacteria“ and „BioFix® Lumi Control solution“ can be stored in a thawed out state in the refrigerator at temperatures of between +2 °C and +8 °C until the expiration date which is printed on them.

The freeze-dried luminescent bacteria should only be removed from the freezer compartment immediately prior to their reactivation. Reactivated luminescent bacteria should be used within 4 hours! An intermediate storage of the reactivated luminescent bacteria should only take place in the refrigerator at temperatures of between +2 °C and +8 °C.

With the advancing storage period of the reactivated luminescent bacteria there can be a reduction of the natural light intensity and a shift of the sensitivity spectrum. The refreezing of reactivated luminescent bacteria cannot be recommended and is not subject to the guarantee obligations of MACHEREY-NAGEL.

Reactivation and dosage of freeze-dried BioFix® Lumi „Single-Shot“ luminescent bacteria:

The reactivation and dosage of freeze-dried BioFix® Lumi „Single-Shot“ luminescent bacteria is always carried out in the same way irrespective of the type of test. Further details to the test procedures can be derived from the instructions for use of the corresponding tests and test systems.

Step 1: Reactivation

1. Remove a deep frozen vial with BioFix® Lumi luminescent bacteria from the deep-freeze compartment and the precooled flask with „BioFix® Lumi Reactivation solution for „Single-Shot“ luminescent bacteria“ from the refrigerator.
2. As quickly as possible add **1,0 mL „BioFix® Lumi Reactivation solution for „Single-Shot“ luminescent bacteria“** („shock thawing“)
3. Mix the vial carefully with reactivated BioFix® Lumi „Single-Shot“ luminous bacteria and label the vial with the word „control“.

Step 2: Dosage

1. Prepare an empty glass cuvette, 50 × 12 mm and label it with the word „sample“. Following this, pipette **0,5 mL of the reactivated luminous bacteria suspension** in the glass cuvette which is labelled with the word „sample“.
2. Leave the „sample“ and the „control“ cuvette to stand for about 10 minutes at + 15 °C.
3. Start the test by measuring the initial luminescence I_0 .
Further details to the test procedure can be derived from the respective instructions for use of the luminometers used.

Analytical quality assurance:

The observance of the MACHEREY-NAGEL internal quality standards which have been implemented in accordance with ISO 9001 is guaranteed with the enclosed test certificate. According to the above mentioned norm the sensitivity of the luminescent bacteria has to be controlled by the end user. The required information such as reference substances of concentrations can be derived from the enclosed test certificate.

pH-value:

According to the DIN EN ISO 11348-3 no correction of the pH is necessary if the pH value of the sample is between 6.0 and 8.5. It may be noted that pH-values of less than 6.0 and greater than 8.5 can lead to pH-related inhibitions of luminescence.

Interferences:

Non-soluble or volatile substances, compounds which are difficult to dissolve, or substances which react with the diluent or the test suspension or which change their state during the test can falsify the test result or make it difficult to reproduce the test results.

Strong colored or turbid samples can cause losses of light due to light adsorption or scattering. In accordance with the procedure described in DIN EN ISO 11348-3, annex A, these effects can be minimized by an appropriate sample treatment (centrifugation or sedimentation) or by using color correction cuvettes (REF 940006).

Oxygen is required for the bioluminescence (> 3 mg/L) and can be limiting in case of samples with a strong oxygen consumption. This can finally result in an inhibition of light emission. If the water sample contains easily biodegradable nutrients a loss of the light intensity which is not caused by contaminants can occur. As the test organism *Vibrio fischeri* is a marine organism the use of sea-water samples can lead to an increase of luminescence masking potential inhibitory effects of the sample (see DIN EN ISO 11348 annex D).

Salt concentrations of the sample higher than 30 g/L NaCl or concentrations of other compounds with similar osmolarity can in conjunction with the salts in the test solutions lead to hyperosmotic effects. To prevent these effects the overall salt concentration should not exceed the osmolarity corresponding to 35 g/L NaCl.

Chlorine containing samples effect the test procedure so that a dechlorination is mandatory, e.g., by addition of a 1 % sodium thiosulfate solution.

Disposal:

Luminescent bacteria and test preparations can be disposed of without difficulty via the sink. Restrictions can arise if the sample contains hazardous or toxic substances which are subject to waste disposal regulations. The user is him/herself responsible for the orderly disposal of such test preparations in accordance with the respectively valid guidelines and regulations.

Literature:

DIN EN ISO 11348-3:2009-05

Determination of the inhibitory effect of water samples upon the light emission of *Vibrio fischeri* (luminescent bacteria test); part 3; procedure with freeze-dried bacteria.

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