

Customer information

Parameter update for compact photometer PF-12^{Plus}

Overview	
Photometer	Compact photometer PF-12 ^{Plus}
Parameter version	(17)
Availability	from 20 th April 2026
Price	free of charge
Link	https://www.mn-net.com/de/nanocolor-device-updates
Required accessories	USB-AB-Connection cable (REF 919687) Computer



Update procedure

NOTICE
Before updating the parameters, it is essential to update the firmware of the PF-12 ^{Plus} to a firmware version $\geq 3.0.1$. Otherwise, the new NANOCOLOR [®] ECO tests will not be displayed correctly and individual methods may no longer be available.

- The parameter file of the PF-12^{Plus} is updated with the parameter update software. An instruction manual on using this software can be found in the download package, which is available via the above mentioned link.
- After a successful parameter update, the following version number will be shown during the device start behind the firmware version in brackets: **(17)**

Modifications

- Implementation of the new NANOCOLOR[®] Permanganate Index 10 (REF 985016)
- Optimization of Method 0-25 NANOCOLOR[®] easily liberated cyanide 04 (REF 985025)

Contact

If you have further questions, please do not hesitate to contact us:

Technical support and customer service	Tel.:	+49 24 21 969-333
	E-Mail:	support@mn-net.com
For orders (international)	E-Mail:	sales@mn-net.com