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Tel.: +1 484 821 0984  
Fax: +1 484 821 1272  
E-mail: sales-us@mn-net.com

## 1. Introduction

The following instruction describes in detail how to update your QUANTOFIX® Relax or URYXXON® Relax easily and quickly. You can update your instrument via Update-Software (3. Relax – Update via Update-Software) or from instrument to instrument (4. Relax – Update from instrument to instrument).

## 2. Special Notes

Please read the following notes carefully before performing an update:

- Do not disconnect the connection between instrument to instrument or computer to instrument during an update. Otherwise, this can cause damage to your instrument.
- Do not switch off the instrument during an update process. Otherwise, this can cause damage to your instrument.
- Please consider that the instrument should only be updated when connected via mains adapter. Do not update the instrument in battery operation.
- The update and subsequent "Load Default" can erase all entries from the memory.
- If you intend an update from instrument to instrument, please consider that this is only possible with the following serial number constellations (please find the serial number on the bottom of the instrument):

Serial number instrument A		Serial number instrument B
URYXXON® Relax	UR2xxxx	UR2xxxx
URYXXON® Relax	URxxxx	URxxxx
QUANTOFIX® Relax	QR2xxxx	QR2xxxx
QUANTOFIX® Relax	QRxxxx	QRxxxx

## 3. Relax – Update via Update-Software

The Update-Software for Relax instruments allows you to update your QUANTOFIX® Relax or URYXXON® Relax easily and quickly.

### Precondition:

Operating system Microsoft® Windows 7, Windows 8 or Windows 10

### Required equipment:

- Instrument: QUANTOFIX® / URYXXON® Relax which needs to be updated
- Update-Software for Relax instruments (Please contact MACHEREY-NAGEL or your local distributor)
- Update-File, e. g. relax.prg (This file is only needed, if no internet connection is available)
- RS232 connection cable (REF 919 772)

### Notes:

The update requires a connection between the computers' and the instruments' RS232 connection. It is not possible to update the instrument via the USB-port. If your computer does not have a natural RS232 port, please make use of a USB-to-RS232 adaptor cable (REF 919 926).

In case of internet availability during the update process, the Update-Software will download the Update-File from the internet automatically. Step number 10 from the instruction below can then be neglected. In case that there is no connection to the internet, you can either use the Update-File stored in the software or select the Update-File manually as described in step number 10 of the instruction.

### Instruction:

1. Open the folder "Update Software" on your computer.
2. Connect the instrument to a power supply.
3. Connect the instrument via the RS232 connection cable to your computer (Please observe information of your operating system, e. g. drivers not found.).
4. Turn on the instrument and tap on the screen several times immediately.
5. Please enter the service code 1234 and confirm your entry with ☒.
6. Go to "Program update" and confirm your entry with ☒.
7. Start the software on the computer by using the "Relax update" icon.
8. In the software screen, please press the binoculars (🔍) for searching and connecting the instrument.
9. Once the instrument has been found, the software displays COM-Port and software version.
10. Optional: The folder icon (📁) becomes active. Please press  and select the Update-File (e. g. relax.prg) in the respective directory.
11. The software confirms that a suitable file has been found and the checkmark (✅) becomes active.
12. Please press ☒ to start the update.
13. Once the update is performed, the instrument automatically restarts and performs a self-check.

www.mn-net.com

MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany

DE / International:

Tel.: +49 24 21 969-0

Fax: +49 24 21 969-199

E-mail: info@mn-net.com

CH:

Tel.: +41 62 388 55 00

Fax: +41 62 388 55 05

E-mail: sales-ch@mn-net.com

FR:

Tel.: +33 388 68 22 68

Fax: +33 388 51 76 88

E-mail: sales-fr@mn-net.com

US:

Tel.: +1 484 821 0984

Fax: +1 484 821 1272

E-mail: sales-us@mn-net.com

14. To finish the procedure, please perform a reset. To do this, switch off the instrument and repeat steps 4 and 5.
15. Go to "Load Default" and confirm your entry with ☒. The screen refreshes and the reset has been performed.
16. Leave the service menu with ☐ X.
17. The update has been performed and the Relax is ready to use with the new firmware version.
18. The new version will be displayed during instrument is started and on the printout.
19. Disconnect the device from the computer and close the Update-Software.

## 4. Relax – Update from instrument to instrument

You can update your QUANTOFIX® Relax or URYXXON® Relax without a computer easily and quickly. You simply need an additional instrument that already contains the desired firmware version.

### Required equipment:

- **Instrument A:** QUANTOFIX® / URYXXON® Relax with new/desired firmware version as a "master instrument"
- **Instrument B:** QUANTOFIX® / URYXXON® Relax which needs to be updated
- RS232 connection cable with accessories (REF 919 774)

### Instruction:

1. Please label **instrument A** and **instrument B** to avoid mistakes during the update process.
2. Please connect **instrument A** and **instrument B** to a power source.
3. Connect **instrument A** to **instrument B** with the RS232 connection cable.
4. Turn on **instrument B** and tap on the screen several times immediately.
5. Please enter the service code 1234 and confirm your entry with ☒.
6. Go to "Program update" and confirm your entry with ☒.
7. Turn on **instrument A** and tap on the screen several times immediately.
8. Please enter the service code 2222 and confirm your entry with ☒.
9. The screen of **instrument B** goes dark, **instrument A** shows the information "connected" and displays the update progress.
10. Once the update is performed, **instrument B** automatically restarts and performs a self-check.
11. To finish the procedure, please perform a reset for **instrument B**. To do this, switch off the **instrument B** and repeat steps 4 and 5.
12. Go to "Load Default" and confirm your entry with ☒. The screen of **instrument B** refreshes and the reset has been performed.
13. Leave the service menu with ☐ X.
14. The update process from **instrument B** has been performed and the Relax is ready to use with the new firmware version.
15. The new version of **instrument B** will be displayed during instrument is started and on the printout.
16. Disconnect both instruments and switch off **instrument A**.

## 5. Contact

For further questions, please do not hesitate to contact us:

<b>Technical Support and Customer Service</b>	Tel.:	+49 24 21 969-332
	E-Mail:	csc@mn-net.com
<b>For orders (international)</b>	Tel.:	0800 2616 000
	E-Mail:	sales-de@mn-net.com

www.mn-net.com

MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany

DE / International:

Tel.: +49 24 21 969-0

Fax: +49 24 21 969-199

E-mail: info@mn-net.com

CH:

Tel.: +41 62 388 55 00

Fax: +41 62 388 55 05

E-mail: sales-ch@mn-net.com

FR:

Tel.: +33 388 68 22 68

Fax: +33 388 51 76 88

E-mail: sales-fr@mn-net.com

US:

Tel.: +1 484 821 0984

Fax: +1 484 821 1272

E-mail: sales-us@mn-net.com