

Customer information

Parameter update for compact photometer PF-12^{Plus}

Overview	
Photometer	Compact photometer PF-12 ^{Plus}
Parameter version	(15)
Availability	from 10 th July 2024
Price	free of charge
Link	https://www.mn-net.com/de/nanocolor-device-updates
Required accessories	USB-AB-Connection cable (REF 919687) Computer



Update procedure

NOTICE

Before updating the parameters, it is essential to update the firmware of the PF-12^{Plus} to a firmware version $\geq 3.0.1$. Otherwise, the new NANOCOLOR[®] ECO tests will not be displayed correctly and individual methods may no longer be available.

- The parameter file of the PF-12^{Plus} is updated with the parameter update software. An instruction manual on using this software can be found in the download package, which is available via the above mentioned link.
- After a successful parameter update, the following version number will be shown during the device start behind the firmware version in brackets: **(15)**

Modifications

- Implementation of the new NANOCOLOR[®] test Boron 3 (REF 985013)
- The test NANOCOLOR[®] Chloride 200 no longer requires a ZERO measurement (REF 985019)
- Optimization of the VISOCOLOR[®] Powder Pillows pH method (REF 936222)
- Bugfixes

Contact

If you have further questions, please do not hesitate to contact us:

Technical support and customer service

Tel.: +49 24 21 969-333

E-Mail: support@mn-net.com

For orders (international)

E-Mail: sales@mn-net.com