MACHEREY-NAGEL

Parameter update software

NANOCOLOR[®] PF-12^{Plus}

- Bedienungsanleitung
- Operating instruction
- Mode d'emploi

Inhalt

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E-mail: sales-de@mn-net.com

1. Introduction

The parameter-updatesoftware for the compact photometer PF-12^{Plus} allows you to update your PF-12^{Plus} easily and quickly with the latest set of parameters.

2. Safety instructions

BE SURE TO READ THE FOLLOWING SAFETY INSTRUCTIONS CAREFULLY BEFORE USING THE SOFTWARE.

Failure to follow these instructions may result in malfunction or damage of the device.

NOTICE

Indicates a situation which, if not avoided, may result in malfunction or damage to the device. Information that clarifies facts in the text and that requires special attention.

3. Precondition

- Operating system Microsoft® Windows 10 / 8 / 7 / Vista
- Instrument PF-12^{Plus} from firmware-version 1.0.6 onwards

4. Required equipment

- Instrument: PF-12^{Plus}
- Parameter-updatesoftware for the PF-12^{Plus}
www.mn-net.com/de/nanocolor-device-updates
- Parameter update file (Included in the download package of the parameter update software. Only needed when the connection to the MN-webserver fails.)
- NANOCOLOR® USB cable A/B (REF 919687). Included in the scope of delivery of the device.

5. Note

- Do not interrupt the connection between the computer and the instrument during the update process. Otherwise, this may cause damage to your instrument.
- Unusable control elements are displayed blurred in the parameter-updatesoftware.

6. Run the software

1. Unzip the folder after the download.
2. Open the folder that includes the software files.
3. Double-click on Method-Update

7. Explanation of the symbols

Symbol	Meaning
	Searches for a connected PF-12 ^{Plus} and establishes the connection.
	Enables the parameter update to be loaded from a local data carrier.
	Starts the parameter update.
	Terminates the program.

8. Perform parameter update

1. Connect the instrument via the USB cable to your computer.
2. Turn on the instrument.
3. Start the software as described in chapter 3.

NOTICE

To establish a successful connection, the device must be switched on and the start process must be completed.

4. Please press  to search for a connected instrument. All instruments which have been found are shown in a drop-down list. Please select the desired instrument on the basis of the serial number (e. g. NPF12P0205) which you can find on the backside of the instrument.

5. Once the instrument has been found, the current *Device Version*, the *Loaded Version* as well as the *Version Source* are shown. The following options are possible in the field Version Source:
 - A The software could not be connected to the MN-webserver. The update will proceed with the updatefile implemented in the software. Attention: This version may be out of date.
 - W The software is connected to the MN-webserver. The update will proceed with the latest updatefile taken from the webserver.
 - F If there is no connection to our webserver, the latest updatefile can be downloaded via www.mn-net.com/de/nanocolor-device-updates and added to the software via .

NOTICE

Do not switch off the device during the update process and do not disconnect computer and device.

6. Please press  to start the parameter update. The progress of the update is indicated by the green status bar.
7. After the update has been performed successfully, the instrument is updated with the new parameter update.

9. Troubleshooting

Error	Source	Solution
Device search without result.	No device was found. The interface of the device is defective. The drivers required for the device are not installed.	Check if the device is connected to the computer and turned on. Check in the device manager whether the necessary drivers are present and installed them manually if necessary. Check if the connection cable is damaged. Check the interface of the device.
No data were found on the server or the server is unreachable.	The software used is out of date. There is no connection to the MACHEREY-NAGEL server.	Check if the latest software version is used. Check the internet connection. Manually load the firmware update file and install it by clicking yes in the error message.

Contact

For further questions, please do not hesitate to contact us:

Technical support and customer service

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