REF 963911

04.23

de

NANOCOLOR®

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en

NANOCOLOR[®]

Cartridges for chloride elimination

Method:

Cartridges for chloride elimination, e.g. in COD or nitrate tests

Contents:

10 cartridges for chloride elimination

Hazard warning:

Cartridges contain silver sulfate 10–30 %.

H318 Causes serious eye damage.

P280sh, P305+351+338, P310 Wear protective gloves/eye protection. IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. Immediately call a POISON CENTER/doctor. For further information ask for a safety data sheet.

Interferences:

Chloride contents up to 2000 mg/L are eliminated by one cartridge pretreatment. For the determination of the chloride content we recommend a preliminary test with QUANTOFIX[®] Chloride (REF 91321).

Note:

The cartridge contents are sensitive to heat and light. Thus, it is recommended to open the cartridge packaging just before use. Avoid contact with the cartridge content. In case of contact with skin, rinse thoroughly with water. Slight color changes of the cartridge content do not interfere with the determination.

Procedure:

Requisite accessories: reaction tubes 16 mm OD (REF 91680), alternatively beaker

Open packaging and take out cartridge.

Dip cartridge into test sample and draw up plunger slowly until about 5 mL of sample solution are inside the cartridge. Shake cartridge well for 30 s. Afterwards, press out sample solution carefully; reject the first drops (ca. 0.5 mL) and fill the remaining contents (clear solution) into an empty reaction tube by slowly pouring the solution down the inner side of the tube.

This pretreated and chloride-free solution can be used for the respective determination.

Disposal:

After use, cartridges can be disposed into waste.

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MACHEREY-NAGEL GmbH & Co. KG · Valencienn Str. 11 · 52355 Düren · Allemagne
France : MACHEREY-NAGEL SAS · 1, rue Gutenberg – BP135 · 67720 Hoerd · France
MACHEREY-NAGEL SAS (Société par Actions Simplifiée) au capital de 186600 €