

User manuals

NucleoSpin® 96 Plant II	740663.2 / .4 / .24
NucleoSpin® 96 Plant II Core Kit	740468.4



Dear valued customer,

Thank you for choosing MACHEREY-NAGEL for your application. We have attached a short protocol for your review.

To obtain the best results we recommend to follow the detailed protocol available online, especially when you are a first time user of this kit.

All important links to the above mentioned products are listed in this leaflet.

Important information regarding product components, specifications, safety instructions, and processing protocols, can all be found on the product website and accessed easily via the QR code.

QR-Code product website



[qr.mn-net.com/qr/\(241\)740663](https://qr.mn-net.com/qr/(241)740663)

Use the following QR code or the link below for direct access to the user manual.

QR-Code user manual



[qr.mn-net.com/qr/\(IFU\)740663](https://qr.mn-net.com/qr/(IFU)740663)

It is strongly recommended to read the detailed protocol section of the user manual if using the kit for the first time. However, experienced users may refer to the protocol at a glance. The protocol at a glance is designed to be used only as a supplemental tool for quick reference while performing the purification procedure.

We are constantly improving our products and we reserve the right to make changes or additions to protocols. Please check for updated revisions for previously downloaded manuals.

This leaflet does not replace the full manual!

Protocol at a glance

NucleoSpin® 96 Plant II – centrifuge processing

Please check the user manual if your centrifuge and setup meets the requirements for centrifuge processing of 96-well plates.

1 Homogenize samples	Up to 100 mg wet or 20 mg lyophilized plant tissue 5,600–6,000 x g, 2 min
2a Cell lysis using Buffer PL1	500 µL PL1 10 µL RNase A Mix 65 °C, 30 min Proceed with step 3
2b Cell lysis using Buffer PL2 and PL3	400 µL PL2 10 µL RNase A Mix 65 °C, 30 min 100 µL PL3 Mix and incubate on ice for 5 min Proceed with step 3
3 Clear lysate by centrifugation	5,600–6,000 x g, 20 min
4 Adjust binding conditions	Mix 450 µL PC with 400 µL cleared lysate
5 Transfer lysate to NucleoSpin® Plant II Binding Plate	
6 Bind DNA to silica membrane of the NucleoSpin® Plant II Binding Plate	5,600–6,000 x g, 2 min
7 Wash and dry silica membrane	400 µL PW1 5,600–6,000 x g, 2 min 700 µL PW2 5,600–6,000 x g, 2 min 700 µL PW2 5,600–6,000 x g, 10 min
8 Elute DNA	100 µL PE (70 °C) (incubate 2 min) 5,600–6,000 x g, 2 min Repeat once

Protocol at a glance

NucleoSpin® 96 Plant II – vacuum processing

1 Homogenize samples	Up to 100 mg wet or 20 mg lyophilized plant tissue 5,600–6,000 x g, 2 min
2a Cell lysis using Buffer PL1	500 µL PL1 10 µL RNase A Mix 65 °C, 30 min Proceed with step 3
2b Cell lysis using Buffer PL2 and PL3	400 µL PL2 Mix 65 °C, 30 min 100 µL PL3 Mix and incubate on ice for 5 min Proceed with step 3
3 Clear lysate by centrifugation	5,000–6,000 x g 20 min
4 Adjust binding conditions	Mix 450 µL PC with 400 µL cleared lysate
5 Transfer lysate to NucleoSpin® Plant II Binding Plate	
6 Bind DNA to silica membrane of the NucleoSpin® Plant II Binding Plate	-0.2 to -0.4 bar* (2 min)
7 Wash and dry silica membrane	400 µL PW1 700 µL PW2 700 µL PW2 -0.4 bar* (1 min each step) Remove MN Wash Plate Dry silica membrane (10 min, maximum vacuum)
8 Elute DNA	100 µL PE (incubate 2 min) -0.4 bar* (2 min) Repeat once

* Reduction of atmospheric pressure

QR-Code Safety Data Sheet

Please find a digital version of the safety data sheets by following the link below:



www.mn-net.com/sds



We strongly recommend to carefully read the detailed protocol section of the product's user manual. If you have any questions about the protocol or product, please contact our Technical Support.

Contact MN

Germany and international

MACHEREY-NAGEL GmbH & Co. KG
Valenciennner Str. 11 · 52355 Düren · Germany
Tel.: +49 24 21 969-0
Toll-free: 0800 26 16 000 (Germany only)
E-mail: info@mn-net.com

Technical support Bioanalysis

Tel.: +49 24 21 969-333
E-mail: support@mn-net.com

USA

MACHEREY-NAGEL Inc.
924 Marcon Blvd. · Suite 102 · Allentown PA, 18109 · USA
Toll-free: 888 321 6224 (MACH)
E-mail: sales-us@mn-net.com

France

MACHEREY-NAGEL SAS
1, rue Gutenberg – BP135 · 67720 Hoerdt Cedex · France
Tel.: +33 388 68 22 68
E-mail: sales-fr@mn-net.com

MACHEREY-NAGEL SAS (Société par Actions Simplifiée) au capital de 186600 €

Siret 379 859 531 00020 · RCS Strasbourg B379859531
N° intracommunautaire FR04 379 859 531

Switzerland

MACHEREY-NAGEL AG
Hirsackerstr. 7 · 4702 Oensingen · Switzerland
Tel.: +41 62 388 55 00
E-mail: sales-ch@mn-net.com