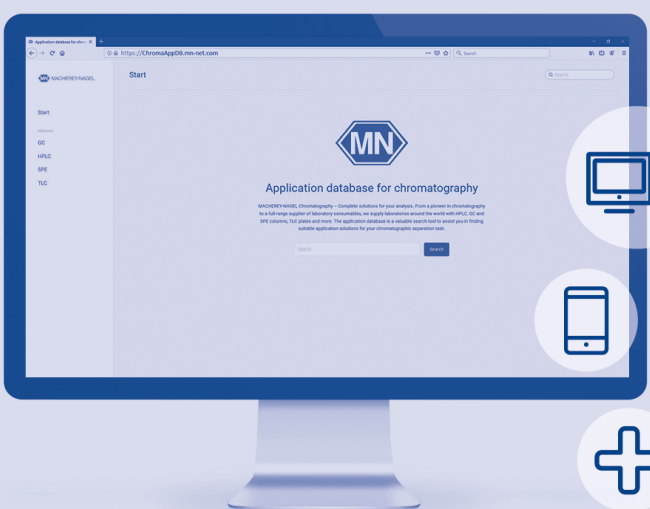
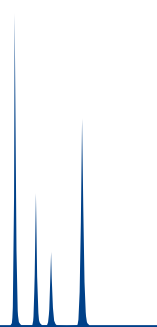
	<b>Column Protection System</b> Vorsäulenhalter – Gebrauchsanleitung	<input type="button" value="de"/>
	<b>Column Protection System</b> Guard column holder – User manual	<input type="button" value="en"/>
	<b>Column Protection System</b> Support de pré-colonne – Manuel	<input type="button" value="fr"/>
	<b>Column Protection System</b> Soporte para columna de protección – Manual de usuario	<input type="button" value="es"/>

Please check the full range of MACHEREY-NAGEL chromatography products: [www.mn-net.com/chromatography](http://www.mn-net.com/chromatography)

 HPLC	 GC	 TLC			
 SPE and Flash	 Syringe filters				
 Vials and caps					
					

... for applicative support please visit our application database with more than 3000 chromatography applications: [ChromaAppDB.mn-net.com](http://ChromaAppDB.mn-net.com) or contact us: [support@mn-net.com](mailto:support@mn-net.com)

**Germany and international:**  
**MACHEREY-NAGEL GmbH & Co. KG**  
 Valenciener Str. 11 · 52355 Düren · Germany  
 Tel.: +49 24 21 969-0  
 info@mn-net.com · [www.mn-net.com](http://www.mn-net.com)

**Switzerland:**  
**MACHEREY-NAGEL AG**  
 Hirsackerstr. 7 · 4702 Oensingen · Switzerland  
 Tel.: 062 388 55 00  
 sales-ch@mn-net.com

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en

REF 718966

The Column Protection System is a universal guard column holder system, suitable for all analytical HPLC columns with 1/16" fittings. It offers:


- ideal protection for your analytical main column → significant increase of column life time
- minimized void volume → suitable also for ultra fast HPLC
- pressure stability up to 1300 bar (18 850 psi)
- visual contamination check → in-time changing of the guard column

MACHEREY-NAGEL offers corresponding EC guard columns for the Column Protection System filled with **NUCLEODUR**<sup>®</sup>, **NUCLEOSHELL**<sup>®</sup>, **NUCLEOSIL**<sup>®</sup> and further HPLC adsorbents. Please check [www.mn-net.com](http://www.mn-net.com) for a detailed description of our modern HPLC phases and a listing of EC guard columns corresponding to your analytical main column.


### Content Column Protection System (REF 718966)

- 1 x Guard column holder
- 2 x Capillaries (ID 0.12 mm)
- 3 x Ferrules (for HPLC columns with particle size > 2 µm)
- 3 x Ferrules (for HPLC columns with particle size < 2 µm)
- 2 x Wrenches (wrench size: 12 and 14 mm)
- 1 x User manual

### Construction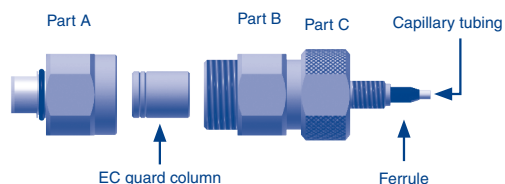


### Guard column selection

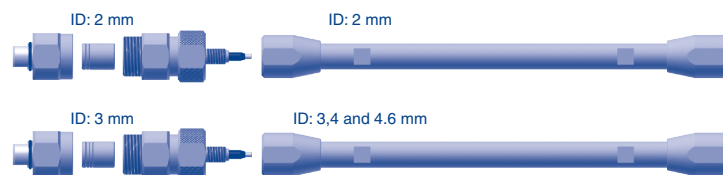
### Installation

#### Step 1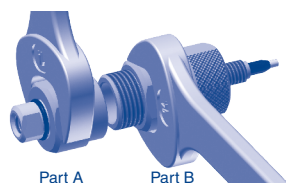


Open the guard column holder (part A and B) by using the wrenches (size 14 mm).

#### Step 2

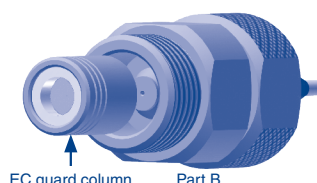
Insert the guard column into the guard column holder part B. We recommend to insert the guard column as shown in the figure.

#### Step 3

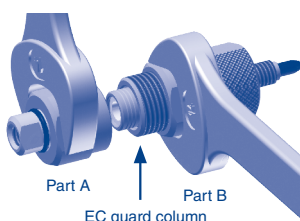Close the guard column holder (part A and B) with the wrenches (size 14 mm).

#### Step 4

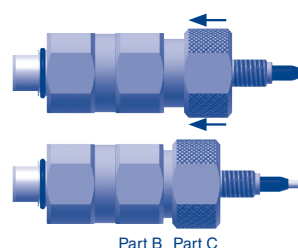Press part C into part B.

#### Step 5


Connect the guard column holder (part C) with the HPLC column by using the wrenches (size 14 and 12 mm).

#### Step 6

Connect the injector capillary with the guard column holder and the HPLC column outlet with the detector capillary.


### EC guard column replacement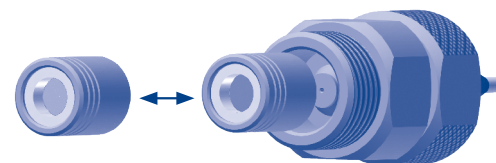


It is advisable to replace the guard column if the silver membrane is contaminated (bright or dark discoloration). If the contaminants are colorless, replace the guard column if the pressure rises or the chromatographic performance decreases.

### Maintenance

#### A) Replace the metal capillary tubing if it is damaged

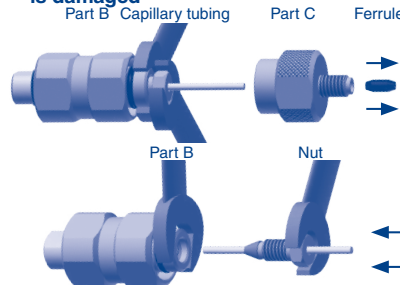#### B) Replace the ferrule in order to prevent leaking
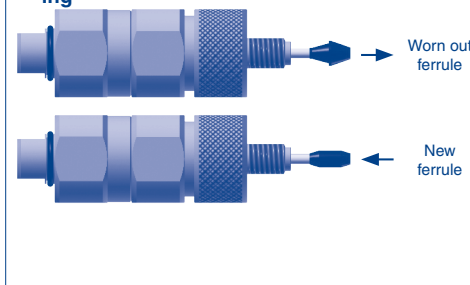
#### C) Replace part C if the thread is worn out


### Troubleshooting

#### Bad chromatographic performance

- Make sure you seated the tubing from the injector all the way down into the entrance of the guard column holder (especially if you use metal capillary tubings).
- Make sure the tubing was cut rectangular.
- Make sure you seated the capillary tubing of the guard column holder all the way down into the entrance of the HPLC column.
- Make sure you wrench tightened the guard column holder units (part A and B) with the wrenches (size: 14 mm) after hand tightening.
- The odds are that the guard column is worn out. Replace the guard column by a new one.

#### Leakage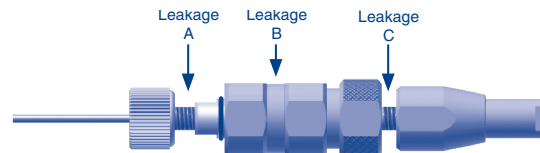

#### Leakage A:

- Make sure that the PEEK fingertight or metal male nut, respectively is not worn out.
- Make sure that the metal capillary has been geared to the inlet of part A of the guard column holder.

#### Leakage B:

- Make sure that you wrench tightened the guard column holder units (part A and B) after hand tightening.
- Make sure the guard column is seated correctly in part B of the guard column holder.

#### Leakage C:

- Make sure that the PEEK ferrule on the metal capillary tubing of the guard column holder is not worn out. If necessary, replace the PEEK ferrule (see maintenance: B).
- Make sure that you seated the capillary tubing of the holder all the way down into the entrance of the HPLC column.
- Make sure that you have fixed the connection between the guard column holder (part C) and the HPLC column with the wrenches

### Replacement parts

- Ferrules for HPLC columns with particle size > 2 µm (pack of 5) (REF 718967)
- Ferrules for HPLC columns with particle size < 2 µm (pack of 5) (REF 718963)
- Replacement connector (part C) including O-ring (pack of 1) (REF 718968)
- Capillary tubings ID 0.12 mm, nuts and metal ferrules (pack of 3) (REF 718969)
- Capillary tubings ID 0.18 mm (for higher flow rates), nuts and metal ferrules (pack of 3) (REF 718971)
- Wrench (size 12 and 14 mm) (pack of 1) (REF 718970)

### UNIVERSAL Guard Columns (for HPLC columns from other manufacturers)

- EC 4/2 UNIVERSAL RP (for HPLC columns with 2 mm ID) (pack of 3) (REF 728777.20)
- EC 4/2 UNIVERSAL RP guard column (for HPLC columns with 2 mm ID), value pack (pack of 9) (REF 728778.20)
- EC 4/3 UNIVERSAL RP (for HPLC columns with 3, 4 or 4.6 mm ID) (pack of 3) (REF 728777.30)
- EC 4/3 UNIVERSAL RP guard column (for main columns with 3, 4 and 4.6 mm ID), value pack (pack of 9) (REF 728778.30)

For **NUCLEODUR**<sup>®</sup>, **NUCLEOSHELL**<sup>®</sup> and **NUCLEOSIL**<sup>®</sup> EC cartridges, please visit [www.mn-net.com](http://www.mn-net.com).

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**MACHEREY-NAGEL SAS**
MACHEREY-NAGEL SAS (Société par Actions Simplifiée)
**MACHEREY-NAGEL Inc.**