

NucleoMag[®] RNA Blood – NucleoMag[®] X32

Protocol details

Application	RNA from stabilized blood
Kit	NucleoMag [®] RNA Blood
REF	744352 (.1 / .4)
Protocol name	NMRNABloodR02a, NMRNABloodR02b
Run time	approx. 78 min



Protocol steps

Procedure
1 Perform lysis according to the user manual NucleoMag [®] RNA Blood.
2 Fill the 96-well Deep-well plates according to the table below.
3 Load the plates on the NucleoMag [®] X32.
4 Insert tip combs on the mounting grooves.*
5 Select the protocol from the instrument menu and start protocol NMRNABloodR02a.
6 After the run (approx. 44 min), dispense 250 µL of Binding Buffer MRB1 to the DNA digestion (column 4 + 10).**
7 Place the plate back into the NucleoMag [®] X32, slide in the tip combs and proceed with protocol NMRNA BloodR02b (approx. 34 min).

Notes: * Please equip all tip combs in order to cover the magnetic rods in used and unused wells.

** To remove the plate the tip combs have to be temporarily removed, they can be reused in their respective positions, but we recommend using a new set of tip combs after placing the plate back into the instrument.

Required consumables and instrumentation provided by user

Product	Specification	REF
NucleoMag [®] X32	Automated nucleic extraction system for MACHEREY-NAGEL's NucleoMag [®] kits enabling parallel processing of up to 32 samples	747020
96 Deep-well Plate	96 deep-well plates for NucleoMag [®] X32 (25 pieces)	744955
Tip Combs	8-well tip combs for NucleoMag [®] X32 (50 pieces)	744960

Find more NucleoMag[®] X32 protocol information and scripts here: <https://www.mn-net.com/de/NucleoMag-X32-Scripts>




Loading table

Position	Reagents	Samples per plate
Column 1 + 7	Sample (820 µL), Proteinase K (15 µL), Binding Buffer MRB1 (140 µL), NucleoMag® B-Beads (15 µL)	Sample 1-8 Sample 9-16
Column 2 + 8	Wash Buffer MRB2 (900 µL)	Sample 1-8 Sample 9-16
Column 3 + 9	Wash Buffer EtOH 80% (900 µL)	Sample 1-8 Sample 9-16
Column 4 + 10	rDNase reaction mixture (300 µL)	Sample 1-8 Sample 9-16
Column 5 + 11	Wash Buffer EtOH 80% (900 µL)	Sample 1-8 Sample 9-16
Column 6 + 12	RNAse-free H ₂ O (100 µL)	Sample 1-8 Sample 9-16

Notes:

Please refer to the image below for a visual representation of the loading scheme

Loading scheme

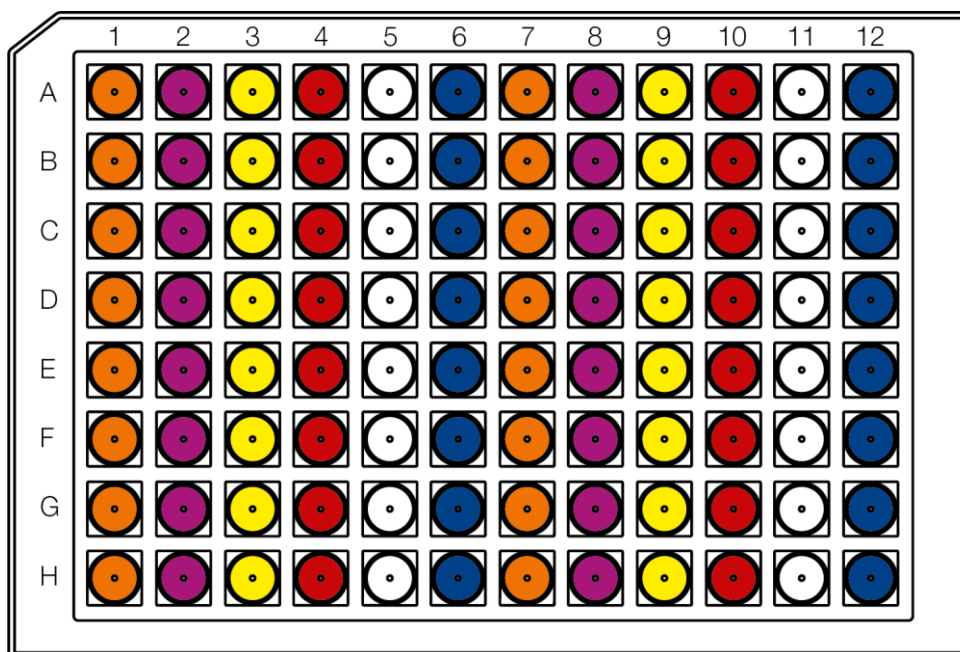

Disclaimer

Information

MACHERY-NAGEL GmbH & Co. KG makes every effort to include accurate and up-to-date information within this publication; however, it is possible that omissions or errors might have occurred. MACHERY-NAGEL GmbH & Co. KG cannot, therefore, make any representations or warranties, expressed or implied, as to the accuracy or completeness of the information provided in this publication. Changes in this publication can be made at any time without notice. For technical details and detailed procedures of the specifications provided in this document please contact your MACHERY-NAGEL representative. This publication may contain reference to applications and products which are not available in all markets. Please check with your local sales representative. All mentioned trademarks are protected by law. All used names and denotations can be brands, trademarks, or registered labels of their respective owner – also if they are not special denotation. To mention products and brands is only a kind of information (i.e., it does not offend against trademarks and brands and can not be seen as a kind of recommendation or assessment). Reference herein to any specific commercial product, process, or service by trade name, trademark, manufacturer, or otherwise, does not necessarily constitute or imply its endorsement, recommendation, or support by MACHERY-NAGEL GmbH & Co. KG. Any views or opinions expressed herein by the authors' do not necessarily state or reflect those of MACHERY-NAGEL. NucleoMag® is a registered trademark of MACHERY-NAGEL GmbH & Co. KG, Düren, Germany.