

en QuantifyGo® – Scan background

Instruction

To ensure the most accurate analysis of your test strip, please follow these simple guidelines when capturing the image with your mobile phone.

A clear and properly framed photograph is essential for the subsequent digital evaluation:

1. Centering:

The test strip must be placed exactly in the center of the black area provided for this purpose so that it does not touch the top edge or the side edges.

2. Orientation:

The test pad(s) must be facing upward and positioned towards the top side of the image.



3. Markers:

All five markers must be fully visible and completely captured within the image frame.

4. Focus:

The image must be sharp and properly focused.

❗ *When printing this file, make sure that the print size is set to 100%. To check whether the size is correct, you can measure the marks.*

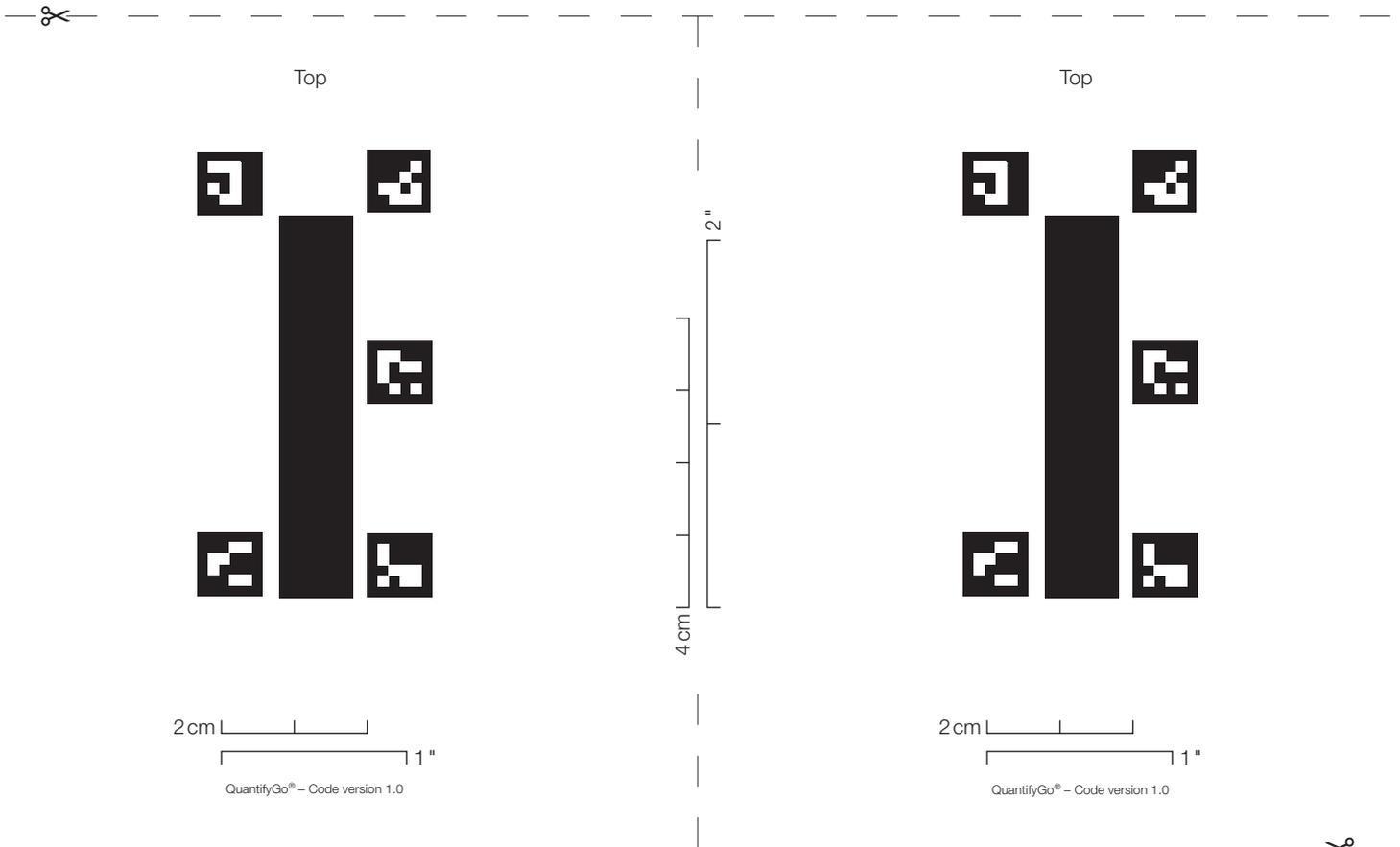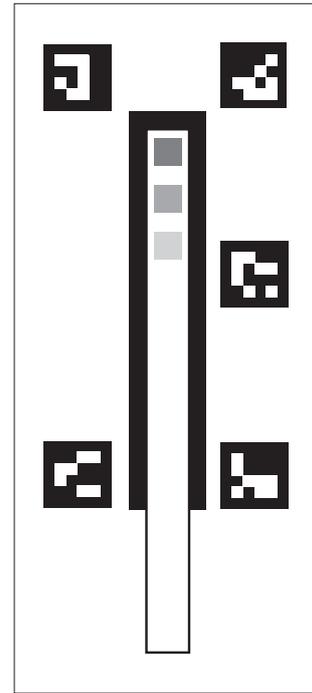

MACHEREY-NAGEL

www.mn-net.com

MACHEREY-NAGEL GmbH & Co. KG · Valencienner Str. 11 · 52355 Düren · Germany

DE +49 24 21 969-0 info@mn-net.com

CH +41 62 388 55 00 sales-ch@mn-net.com

FR +33 388 68 22 68 sales-fr@mn-net.com

US +1 888 321 62 24 sales-us@mn-net.com

